



碧云天生物技术/Beyotime Biotechnology  
订货热线: 400-1683301或800-8283301  
订货e-mail: order@beyotime.com  
技术咨询: info@beyotime.com  
网址: http://www.beyotime.com

## Hoechst 33342

产品编号	产品名称	包装
C1022	Hoechst 33342	10mg

### 产品简介:

- Hoechst 33342, 也称bisBenzimide H 33342或HOE 33342。分子式为C<sub>27</sub>H<sub>28</sub>N<sub>6</sub>O·3HCl·3H<sub>2</sub>O, 分子量为615.99, CAS Number 23491-52-3。
- Hoechst 33342是一种可以穿透细胞膜的蓝色荧光染料, 对细胞的毒性较低。
- Hoechst 33342染色常用于细胞凋亡检测, 染色后用荧光显微镜观察或流式细胞仪检测。Hoechst 33342也常用于普通的细胞核染色, 或常规的DNA染色。
- Hoechst 33342的最大激发波长为346nm, 最大发射波长为460nm; Hoechst 33342和双链DNA结合后, 最大激发波长为350nm, 最大发射波长为461nm。
- Hoechst 33342溶于水, 溶解度可达20mg/ml。
- 用于细胞核染色时, 推荐的Hoechst 33342工作浓度为0.5-10μg/ml。

### 包装清单:

产品编号	产品名称	包装
C1022	Hoechst 33342	10mg
—	说明书	1份

### 保存条件:

-20°C避光保存, 一年有效。

### 注意事项:

- 荧光染料都存在淬灭的问题, 建议染色后尽量当天完成检测。
- 为减缓荧光淬灭可以使用抗荧光淬灭封片液。抗荧光淬灭封片液(P0126)可以向碧云天订购。
- Hoechst 33342对人体有害, 操作时请小心, 并注意有效防护以避免直接接触人体或吸入体内。
- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

### 相关产品:

产品编号	产品名称	包装
C1002	DAPI	5mg/ml×0.2ml
C1005	DAPI染色液	10ml
C1006	DAPI染色液	50ml
C1011	Hoechst 33258	10mg
C1017	Hoechst 33258染色液	10ml
C1018	Hoechst 33258染色液	50ml
C1022	Hoechst 33342	10mg
C1025	Hoechst 33342染色液	10ml
C1026	Hoechst 33342染色液	50ml
C1027	Hoechst 33342活细胞染色液(100X)	0.1ml
C1028	Hoechst 33342活细胞染色液(100X)	0.5ml
C1029	Hoechst 33342活细胞染色液(100X)	3ml

### 使用本产品的文献:

1. Li HG, Liao AH, Ding XF, Zhou H, Xiong CL. The expression and significance of CATSPER1 in human testis and ejaculated spermatozoa. Asian J Androl. 2006 May;8(3):301-6.
2. Yao G, Ling L, Luan J, Ye D, Zhu P. Nonylphenol induces apoptosis of Jurkat cells by a caspase-8 dependent mechanism. Int Immunopharmacol. 2007 Apr;7(4):444-53.
3. Wang Q, Wang Y, Liang C, Song J, Chen X. Identification of a hydrophobic domain of HA2 essential to morphogenesis of Helicoverpa armigera nucleopolyhedrovirus. J Virol. 2008 Apr;82(8):4072-81.

4. Badraoui R, Blouin S, Moreau MF, Gallois Y, Rebai T, Sahnoun Z, Baslé M, Chappard D. Effect of alpha tocopherol acetate in Walker 256/B cells-induced oxidative damage in a rat model of breast cancer skeletal metastases. *Chem Biol Interact.* 2009;182(2-3):98-105.
5. Han M, He CX, Fang QL, Yang XC, Diao YY, Xu DH, He QJ, Hu YZ, Liang WQ, Yang B, Gao JQ. A novel camptothecin derivative incorporated in nano-carrier induced distinguished improvement in solubility, stability and anti-tumor activity both in vitro and in vivo. *Pharm Res.* 2009 Apr;26(4):926-35.
6. Chen Z, Li Z, Peng G, Chen X, Yin W, Kotlikoff MI, Yuan ZQ, Ji G. Extracellular ATP-induced nuclear Ca<sup>2+</sup> transient is mediated by inositol 1,4,5-trisphosphate receptors in mouse pancreatic beta-cells. *Biochem Biophys Res Commun.* 2009 May 1;382(2):381-4.
7. Liu Y, Sheng Z, Liu H, Wen D, He Q, Wang S, Shao W, Jiang RJ, An S, Sun Y, Bendena WG, Wang J, Gilbert LI, Wilson TG, Song Q, Li S. Juvenile hormone counteracts the bHLH-PAS transcription factors MET and GCE to prevent caspase-dependent programmed cell death in Drosophila. *Development.* 2009 Jun;136(12):2015-25.
8. Ma K, Hu M, Qi Y, Qiu L, Jin Y, Yu J, Li B. Structure-transfection activity relationships with glucocorticoid-polyethyl-enimine conjugate nuclear gene delivery systems. *Biomaterials.* 2009 Aug;30(22):3780-9.
9. Zeng X, Nan F, Liang C, Song J, Wang Q, Vlak JM, Chen X. Functional analysis of the *Autographa californica* nucleopolyhedrovirus IAP1 and IAP2. *Sci China C Life Sci.* 2009 Aug;52(8):761-70.
10. Ma K, Hu MX, Qi Y, Zou JH, Qiu LY, Jin Y, Ying XY, Sun HY. PAMAM-triamcinolone acetonide conjugate as a nucleus-targeting gene carrier for enhanced transfer activity. *Biomaterials.* 2009 Oct;30(30):6109-18.
11. Gong Y, Wu J, Huang Y, Shen S, Han X. Nonylphenol induces apoptosis in rat testicular Sertoli cells via endoplasmic reticulum stress. *Toxicol Lett.* 2009;186(2):84-95.
12. Liu Y, Huang H, Lin J, Zhang Q, Tan J, Ren J. Suppressive effects of genomic imprinted gene PEG10 on hydrogen peroxide-induced apoptosis in L02 cells. *J Huazhong Univ Sci Technolog Med Sci.* 2009;29(6):705-9.
13. Badraoui R, Blouin S, Moreau MF, Gallois Y, Rebai T, Sahnoun Z, Baslé M, Chappard D. Effect of alpha tocopherol acetate in Walker 256/B cells-induced oxidative damage in a rat model of breast cancer skeletal metastases. *Chem Biol Interact.* 2009;182(2-3):98-105.
14. Chen YJ, Huang XB, Li ZX, Yin LL, Chen WQ, Li L. Tenuigenin protects cultured hippocampal neurons against methylglyoxal-induced neurotoxicity. *Eur J Pharmacol.* 2010;645(1-3):1-8.
15. Zhang L, Li HZ, Gong X, Luo FL, Wang B, Hu N, Wang CD, Zhang Z, Wan JY. Protective effects of Asiaticoside on acute liver injury induced by lipopolysaccharide/D-galactosamine in mice. *Phytomedicine.* 2010;17(10):811-9.
16. Gao X, Zhang X, Yang X. Morphological apoptotic characteristics of the post-meiotic micronuclei in *Paramecium caudatum*. *Eur J Protistol.* 2010;46(3):243-50.
17. Jiang X, Zhang J, Xia M, Qiu W, Wang H, Zhao D, Wang Y. Role of activating transcription factor 3 (ATF3) in sublytic C5b-9-induced glomerular mesangial cell apoptosis. *Cell Mol Immunol.* 2010;7(2):143-51.
18. Wu H, Lin Y, Shi Y, Qian W, Tian Z, Yu Y, Huo K. SIAH-1 interacts with mammalian polyhomeotic homologues HPH2 and affects its stability via the ubiquitin-proteasome pathway. *Biochem Biophys Res Commun.* 2010;397(3):391-6.
19. Wang CJ, Hu CP, Xu KP, Yuan Q, Li FS, Zou H, Tan GS, Li YJ. Protective effect of selaginellin on glutamate-induced cytotoxicity and apoptosis in differentiated PC12 cells. *Naunyn Schmiedebergs Arch Pharmacol.* 2010;381(1):73-81.
20. Guo XF, Zhu XF, Shang Y, Zhang SH, Zhen YS. A bispecific enediyne-energized fusion protein containing ligand-based and antibody-based oligopeptides against epidermal growth factor receptor and human epidermal growth factor receptor 2 shows potent antitumor activity. *Clin Cancer Res.* 2010 Apr 1;16(7):2085-94.
21. Peng ZG, Fan B, Du NN, Wang YP, Gao LM, Li YH, Li YH, Liu F, You XF, Han YX, Zhao ZY, Cen S, Li JR, Song DQ, Jiang JD. Small molecular compounds that inhibit hepatitis C virus replication through destabilizing heatshock cognate 70 messenger RNA. *Hepatology.* 2010 Sep;52(3):845-53.
22. Yang M, Zhang M, Chen J, Mukherjee R, Zhang L, Lin S, Gu Y. Angiopoietin-1 inhibits mouse glomerular endothelial cell senescence via Tie2 receptor-modulated ERK1/2 signaling. *Am J Nephrol.* 2010; 31(6): 490-500.
23. Zhang YP, Wan P, Wang HQ, Zhao H, Xu YX, Yang R, Zhu CQ. Effect of neuronal excitotoxicity on Munc18-1 distribution in nuclei of rat hippocampal neuron and primary cultured neuron. *Neurosci Bull.* 2011 Jun;27(3):163-72.
24. Wang Y, Hong C, Zhou C, Xu D, Qu HB, Cheng Y. Screening Antitumor Compounds Psoralen and Isopsonalen from *Psoralea corylifolia* L. Seeds. *Evid Based Complement Alternat Med.* 2011;2011:363052.
25. Peng ZG, Zhao ZY, Li YP, Wang YP, Hao LH, Fan B, Li YH, Wang YM, Shan YQ, Han YX, Zhu YP, Li JR, You XF, Li ZR, Jiang JD. Host apolipoprotein B messenger RNA-editing enzyme catalytic polypeptide-like 3G is an innate defense factor and drug target against hepatitis C virus. *Hepatology.* 2011 Apr;53(4):1080-9.
26. Hu XL, Cheng X, Cai L, Tan GH, Xu L, Feng XY, Lu TJ, Xiong H, Fei J, Xiong ZQ. Conditional deletion of NRSF in forebrain neurons accelerates epileptogenesis in the kindling model. *Cereb Cortex.* 2011 Sep;21(9):2158-65.
27. He XN, Su F, Lou ZZ, Jia WZ, Song YL, Chang HY, Wu YH, Lan J, He XY, Zhang Y. Ipr1 gene mediates RAW 264.7 macrophage cell line resistance to *Mycobacterium bovis*. *Scand J Immunol.* 2011 Nov;74(5):438-44.
28. Wu H, Shi Y, Lin Y, Qian W, Yu Y, Huo K. Eukaryotic translation elongation factor 1 delta inhibits the ubiquitin ligase activity of SIAH-1. *Mol Cell Biochem.* 2011 Nov;357(1-2):209-15.
29. Yang H, Taylor HS, Lei C, Cheng C, Zhang W. Hormonal regulation of galectin 3 in trophoblasts and its effects on endometrium. *Reprod Sci.* 2011 Nov;18(11):1118-27.
30. Wang H, Xie X, Zhang F, Zhou Q, Tao Q, Zou Y, Chen C, Zhou C, Yu S. Evaluation of cholesterol depletion as a marker of nephrotoxicity in vitro for novel β-cyclodextrin derivatives. *Food Chem Toxicol.* 2011 Jun;49(6):1387-93.
31. Zhou J, Lin J, Zhou C, Deng X, Xia B. Cytotoxicity of red fluorescent protein DsRed is associated with the suppression of Bcl-xL translation. *FEBS Lett.* 2011 Mar 9;585(5):821-7.
32. Liu Q, Kou JP, Yu BY. Ginsenoside Rg1 protects against hydrogen peroxide-induced cell death in PC12 cells via inhibiting NF-κB activation. *Neurochem Int.* 2011 Jan;58(1):119-25.
33. Kong FL, Cheng W, Chen J, Liang Y. d-Ribose glycates β(2)-microglobulin to form aggregates with high cytotoxicity through a ROS-mediated pathway. *Chem Biol Interact.* 2011 Oct 15;194(1):69-78.
34. Ji SZ, Xiao SC, Luo PF, Huang GF, Wang GY, Zhu SH, Wu MJ, Xia ZF. An epidermal stem cells niche microenvironment created by engineered human amniotic membrane. *Biomaterials.* 2011 Nov;32(31):7801-11.
35. Zhang X, Wang J, Zhou Q, Xu Y, Pu S, Wu J, Xue Y, Tian Y, Lu J, Jiang W, Du D. Brain-derived neurotrophic factor-activated astrocytes produce mechanical allodynia in neuropathic pain. *Neuroscience.* 2011 Dec 29;199:452-60.
36. Li Z, Mi L, Wang PN, Chen JY. Study on the visible-light-induced photokilling effect of nitrogen-doped TiO<sub>2</sub> nanoparticles on cancer cells. *Nanoscale Res Lett.* 2011 Apr 21;6(1):356.
37. Liu SL, Zhang ZL, Sun EZ, Peng J, Xie M, Tian ZQ, Lin Y, Pang DW. Visualizing the endocytic and exocytic processes of wheat germ agglutinin by quantum dot-based single-particle tracking. *Biomaterials.* 2011 Oct;32(30):7616-24.
38. Bai D, Gao Q, Li C, Ge L, Gao Y, Wang H. A conserved TGFβ1/HuR feedback circuit regulates the fibrogenic response in fibroblasts. *Cell Signal.* 2012 Jul;24(7):1426-32.
39. Li B, Xu H, Li Z, Yao M, Xie M, Shen H, Shen S, Wang X, Jin Y. Bypassing multidrug resistance in human breast cancer cells with lipid/polymer particle assemblies. *Int J Nanomedicine.* 2012;7:187-97.
40. Wang S, Tan Y, Lei W, Wang Y, Zhou X, Jia X, Zhang K, Chu L, Liu X, Qian W. Complete eradication of xenograft hepatoma by oncolytic adenovirus ZD55 harboring TRAIL-IETD-Smac gene with broad antitumor effect. *Hum Gene Ther.* 2012 Sep;23(9):992-1002.
41. Zhang J, Zhou F, Wu X, Zhang X, Chen Y, Zha BS, Niu F, Lu M, Hao G, Sun Y, Sun J, Peng Y, Wang G. Cellular pharmacokinetic mechanisms of adriamycin resistance and its modulation by 20(S)-ginsenoside Rh2 in MCF-7/Adr cells. *Br J Pharmacol.* 2012 Jan;165(1):120-34.
42. Zhang KJ, Qian J, Wang SB, Yang Y. Targeting Gene-Viro-Therapy with AFP driving Apoptin gene shows potent antitumor effect in hepatocarcinoma. *J Biomed Sci.* 2012 Feb 9;19:20.

43. He G, Lei W, Wang S, Xiao R, Guo K, Xia Y, Zhou X, Zhang K, Liu X, Wang Y. Overexpression of tumor suppressor TSLC1 by a survivin-regulated oncolytic adenovirus significantly inhibits hepatocellular carcinoma growth. *J Cancer Res Clin Oncol.* 2012 Apr;138(4):657-70.
44. Xiong XY, Guo L, Gong YC, Li ZL, Li YP, Liu ZY, Zhou M. In vitro & in vivo targeting behaviors of biotinylated Pluronic F127/poly(lactic acid) nanoparticles through biotin-avidin interaction. *Eur J Pharm Sci.* 2012 Aug 15;46(5):537-44.
45. Yin QQ, Dong CF, Dong SQ, Dong XL, Hong Y, Hou XY, Luo DZ, Pei JJ, Liu XP. AGEs Induce Cell Death via Oxidative and Endoplasmic Reticulum Stresses in Both Human SH-SY5Y Neuroblastoma Cells and Rat Cortical Neurons. *Cell Mol Neurobiol.* 2012 Nov;32(8):1299-309.
46. Li C, Nie H, Wang M, Su L, Li J, Yu B, Wei M, Ju J, Yu Y, Yan M, Gu Q, Zhu Z, Liu B. MicroRNA-409-3p regulates cell proliferation and apoptosis by targeting PHF10 in gastric cancer. *Cancer Lett.* 2012 Jul 28;320(2):189-97.
47. Guo HF, Li ZS, Dong SW, Chen WJ, Deng L, Wang YF, Ying DJ. Piezoelectric PU/PVDF electrospun scaffolds for wound healing applications. *Colloids Surf B Biointerfaces.* 2012 Aug 1;96:29-36.
48. Zhou L, An XF, Teng SC, Liu JS, Shang WB, Zhang AH, Yuan YG, Yu JY. Pretreatment with the total flavone glycosides of *Flos Abelmoschus manihot* and hyperoside prevents glomerular podocyte apoptosis in streptozotocin-induced diabetic nephropathy. *J Med Food.* 2012 May;15(5):461-8.
49. Yin T, Wang P, Zheng R, Zheng B, Cheng D, Zhang X, Shuai X. Nanobubbles for enhanced ultrasound imaging of tumors. *Int J Nanomedicine.* 2012;7:895-904.
50. Zheng J, Chen Z, Yin W, Miao L, Zhou Z, Ji G. Ryanodine receptors are involved in nuclear calcium oscillation in primary pancreatic  $\beta$ -cells. *Biochem Biophys Res Commun.* 2012 Jun 29;423(2):207-11.
51. Tu Z, Li H, Ma Y, Tang B, Tian J, Akers W, Achilefu S, Gu Y. The enhanced antiproliferative response to combined treatment of trichostatin A with raloxifene in MCF-7 breast cancer cells and its relevance to estrogen receptor  $\beta$  expression. *Mol Cell Biochem.* 2012 Jul;366(1-2):111-22.
52. Zheng X, Tian J, Weng L, Wu L, Jin Q, Zhao J, Wang L. Cytotoxicity of cadmium-containing quantum dots based on a study using a microfluidic chip. *Nanotechnology.* 2012 Feb 10;23(5):055102.
53. Tu Z, Ma Y, Tian J, Li H, Akers W, Achilefu S, Gu Y. Estrogen receptor  $\beta$  potentiates the antiproliferative effect of raloxifene and affects the cell migration and invasion in HCT-116 colon cancer cells. *J Cancer Res Clin Oncol.* 2012 Jul;138(7):1091-103.
54. Li S, Zhang Z, Xue J, Liu A, Zhang H. Cold-inducible RNA binding protein inhibits  $H_2O_2$ -induced apoptosis in rat cortical neurons. *Brain Res.* 2012 Mar 2;1441:47-52.
55. Cheng J, Tang W, Su Z, Guo J, Tong L, Wei Q. Calcineurin subunit B promotes TNF-alpha-induced apoptosis by binding to mitochondria and causing mitochondrial  $Ca^{2+}$  overload. *Cancer Lett.* 2012 Aug 28;321(2):169-78.
56. Zhang G, Lan X, Yen TC, Chen Q, Pei Z, Qin C, Zhang Y. Therapeutic gene expression in transduced mesenchymal stem cells can be monitored using a reporter gene. *Nucl Med Biol.* 2012 Nov;39(8):1243-50.
57. Chen LP, Li ZF, Ping M, Li R, Liu J, Xie XH, Song XJ, Guo L. Regulation of Olig2 during astroglial differentiation in the subventricular zone of a cuprizone-induced demyelination mouse model. *Neuroscience.* 2012 Sep 27;221:96-107.
58. Yu J, Xie X, Zheng M, Yu L, Zhang L, Zhao J, Jiang D, Che X. Fabrication and characterization of nuclear localization signal-conjugated glycol chitosan micelles for improving the nuclear delivery of doxorubicin. *Int J Nanomedicine.* 2012;7:5079-90.
59. Li JZ, Yu SY, Wu JH, Shao QR, Dong XM. Paeoniflorin protects myocardial cell from doxorubicin-induced apoptosis through inhibition of NADPH oxidase. *Can J Physiol Pharmacol.* 2012;90(12):1569-75.
60. Zhang KJ, Zhang J, Wu YM, Qian J, Liu XJ, Yan LC, Zhou XM, Xiao RJ, Wang YG, Cao X, Wei N, Liu XR, Tang B, Jiao XY, Chen K, Liu XY. Complete eradication of hepatomas using an oncolytic adenovirus containing AFP promoter controlling E1A and an E1B deletion to drive IL-24 expression. *Cancer Gene Ther.* 2012 Sep;19(9):619-29.
61. Du J, Tang B, Wang J, Sui H, Jin X, Wang L, Wang Z. Antiproliferative effect of alpinetin in BxPC-3 pancreatic cancer cells. *Int J Mol Med.* 2012 Apr;29(4):607-12.
62. Yao Q, Zhang AM, Ma H, Lin S, Wang XX, Sun JG, Chen ZT. Novel molecular beacons to monitor microRNAs in non-small-cell lung cancer. *Mol Cell Probes.* 2012 Oct;26(5):182-7.
63. Zhang B, Wang Y, Pang X. Enhanced radiosensitivity of EC109 cells by inhibition of HDAC1 expression. *Med Oncol.* 2012 Mar;29(1):340-8.
64. Xie X, Sun A, Zhu W, Huang Z, Hu X, Jia J, Zou Y, Ge J. Transplantation of mesenchymal stem cells preconditioned with hydrogen sulfide enhances repair of myocardial infarction in rats. *Tohoku J Exp Med.* 2012;226(1):29-36.
65. Shen HJ, Shi H, Ma K, Xie M, Tang LL, Shen S, Li B, Wang XS, Jin Y. Polyelectrolyte capsules packaging BSA gels for pH-controlled drug loading and release and their antitumor activity. *Acta Biomater.* 2013 Apr;9(4):6123-33.
66. Peng Y, Zhang J, Xu H, He J, Ying X, Wang Y. Neuroprotective effect of protease-activated receptor-2 in the hypoxia-induced apoptosis of rat RGC-5 cells. *J Mol Neurosci.* 2013 May;50(1):98-108.
67. Li J, Cheng Y, Zhang X, Zheng L, Han Z, Li P, Xiao Y, Zhang Q, Wang F. The in vivo immunomodulatory and synergistic anti-tumor activity of thymosin  $\alpha 1$ -thymopentin fusion peptide and its binding to TLR2. *Cancer Lett.* 2013 Sep 1;337(2):237-47.
68. Zhu W, Shi XL, Xiao JQ, Gu GX, Ding YT, Ma ZL. Effects of xenogeneic adipose-derived stem cell transplantation on acute-on-chronic liver failure. *Hepatobiliary Pancreat Dis Int.* 2013;12(1):60-7.
69. Shen H, Shi H, Xie M, Ma K, Li B, Shen S, Wang X, Jin Y. Biodegradable chitosan/alginate BSA-gel-capsules for pH-controlled loading and release of doxorubicin and treatment of pulmonary melanoma. *J Mater Chem B.* 2013 May;32:3906-17.
70. Lei W, Liu HB, Wang SB, Zhou XM, Zheng SD, Guo KN, Ma BY, Xia YL, Tan WS, Liu XY, Wang YG. Tumor suppressor in lung cancer-1 (TSLC1) mediated by dual-regulated oncolytic adenovirus exerts specific antitumor actions in a mouse model. *Acta Pharmacol Sin.* 2013 Apr;34(4):531-40.
71. Zheng L, Li Y, Cui Y, Yin H, Liu T, Yu G, Lv F, Yang J. Generation of an effective anti-lung cancer vaccine by DTPP-mediated photodynamic therapy and mechanistic studies. *Lasers Med Sci.* 2013 Sep;28(5):1383-92.
72. Han F, Xu Y, Jiang D, Qin Y, Chen H. A BODIPY-derived fluorescent probe for cellular pH measurements. *Anal Biochem.* 2013 Apr 15;435(2):106-13.
73. Liu Y, He G, Wang Y, Guan X, Pang X, Zhang B. MCM-2 is a therapeutic target of Trichostatin A in colon cancer cells. *Toxicol Lett.* 2013 Jul 31;221(1):23-30.
74. Zhou HL, Zheng YJ, Cheng XZ, Lv YS, Gao R, Mao HP, Chen Q. Intercellular Transfer of P-Glycoprotein from the Drug Resistant Human Bladder Cancer Cell Line BIU-87 Does Not Require Cell-to-Cell Contact. *J Urol.* 2013 Sep;190(3):1069-75.
75. Zhu K, Guo C, Xia Y, Lai H, Yang W, Wang Y, Song D, Wang C. Transplantation of novel vascular endothelial growth factor gene delivery system manipulated skeletal myoblasts promote myocardial repair. *Int J Cardiol.* 2013 Oct 3;168(3):2622-31.
76. Wang J, Chen Y, Zhang W, Zheng G, Meng S, Che H, Ke T, Yang J, Chen J, Luo W. Akt Activation Protects Liver Cells from Apoptosis in Rats during Acute Cold Exposure. *Int J Biol Sci.* 2013 May 25;9(5):509-17.
77. Chen Q, Sun J. Implication of Oxidative Stress in Dose-Dependent Toxicity of Silica Nanoparticles in Hepatocytes. *Advanced Materials Research.* 2013;647:42-5.
78. Xu C, Liu C, Huang W, Tu S, Wan F. Effect of Mst1 overexpression on the growth of human hepatocellular carcinoma HepG2 cells and the sensitivity to cisplatin in vitro. *Acta Biochim Biophys Sin (Shanghai).* 2013 Apr;45(4):268-79.
79. Yin T, Wang P, Li J, Zheng R, Zheng B, Cheng D, Li R, Lai J, Shuai X. Ultrasound-sensitive siRNA-loaded nanobubbles formed by hetero-assembly of polymeric micelles and liposomes and their therapeutic effect in gliomas. *Biomaterials.* 2013 Jun;34(18):4532-43.
80. Wang B, Yu T, Dong X, Zhang Z, Song L, Xu Y, Zhang XH. Edwardsiella tarda invasion of fish cell lines and the activation of divergent cell death pathways. *Vet Microbiol.* 2013 May 3;163(3-4):282-9.
81. Wang JJ, Liu YL, Sun YC, Ge W, Wang YY, Dyce PW, Hou R, Shen W. Basic Fibroblast Growth Factor Stimulates the Proliferation of Bone Marrow Mesenchymal Stem Cells in Giant Panda (*Ailuropoda melanoleuca*). *PLoS One.* 2015 Sep 16;10(9):e0137712.
82. Hou L, Li Y, Song H, Zhang Z, Sun Y, Zhang X, Wu K. Protective Macroautophagy Is Involved in Vitamin E Succinate Effects on Human Gastric Carcinoma Cell Line SGC-7901 by Inhibiting mTOR Axis

- Phosphorylation. PLoS One. 2015 Jul 13;10(7):e0132829.
83. Wu X, Zhao B, Cheng Y, Yang Y, Huang C, Meng X, Wu B, Zhang L, Lv X, Li J. Melittin induces PTCH1 expression by down-regulating MeCP2 in human hepatocellular carcinoma SMMC-7721 cells. *Toxicol Appl Pharmacol*. 2015 Oct 1;288(1):74-83.
  84. Gao S, Yu Y, Ma ZY, Sun H, Zhang YL, Wang XT, Wang C, Fan WM, Zheng QY, Ma CL. NMDAR-Mediated Hippocampal Neuronal Death is Exacerbated by Activities of ASIC1a. *Neurotox Res*. 2015 Aug; 28(2): 122-37.
  85. Wang X, Ma Z, Fu Z, Gao S, Yang L, Jin Y, Sun H, Wang C, Fan W, Chen L, Zheng QY, Bi G, Ma CL. Hydroxysafflor Yellow A Protects Neurons From Excitotoxic Death through Inhibition of NMDARs. *ASN Neuro*. 2016 Apr 10;8(2). pii: 1759091416642345.
  86. Yang F, Yan G, Li Y, Han Z, Zhang L, Chen S, Feng C, Huang Q, Ding F, Yu Y, Bi C, Cai B, Yang L. Astragalus Polysaccharide Attenuated Iron Overload-Induced Dysfunction of Mesenchymal Stem Cells via Suppressing Mitochondrial ROS. *Cell Physiol Biochem*. 2016; 39(4): 1369-79.
  87. Li L, Liu JC, Lai FN, Liu HQ, Zhang XF, Dyce PW, Shen W, Chen H. Di (2-ethylhexyl) Phthalate Exposure Impairs Growth of Antral Follicle in Mice. *PLoS One*. 2016 Feb 4;11(2):e0148350.
  88. Wang Y, Qiu Y, Yin S, Zhang L, Shi K, Gao H, Zhang Z, He Q. A functional nanocarrier that copenetrates extracellular matrix and multiple layers of tumor cells for sequential and deep tumor autophagy inhibitor and chemotherapeutic delivery. *Autophagy*. 2016 Dec 2:1-12.
  89. Liu Y, Zhi D, Li M, Liu D, Wang X, Wu Z, Zhang Z, Fei D, Li Y, Zhu H, Xie Q, Yang H, Li H. Shengmai Formula suppressed over-activated Ras/MAPK pathway in *C. elegans* by opening mitochondrial permeability transition pore via regulating cyclophilin D. *Sci Rep*. 2016 Dec 16; 6: 38934.
  90. Liu Y, Zhi D, Li M, Liu D, Wang X, Wu Z, Zhang Z, Fei D, Li Y, Zhu H, Xie Q, Yang H, Li H. Shengmai Formula suppressed over-activated Ras/MAPK pathway in *C. elegans* by opening mitochondrial permeability transition pore via regulating cyclophilin D. *Sci REP-UK*. 2016 Dec 16;6:38934
  91. You F, Li Q, Jin G, Zheng Y, Chen J, Yang H. Genistein protects against A  $\beta$  25-35 induced apoptosis of PC12 cells through JNK signaling and modulation of Bcl-2 family messengers. *BMC Neurosci*. 2017 Jan 12;18(1):12
  92. Wang Z, Liao K, Zuo W, Liu X, Qiu Z, Gong Z, Liu C, Zeng Q, Qian Y, Jiang L, Bu Y, Hong S, Hu G. Depletion of NFBD1/MDC1 Induces Apoptosis in Nasopharyngeal Carcinoma Cells Through the p53-ROS-Mitochondrial Pathway. *Oncol Res*. 2017 Jan 2;25(1):123-136
  93. Hu Y, Xiong LL, Zhang P, Wang TH. Microarray expression profiles of genes in lung tissues of rats subjected to focal cerebral ischemia-induced lung injury following bone marrow-derived mesenchymal stem cell transplantation. *Int J Mol Med*. 2017 Jan;39(1):57-70
  94. Zhang Y, Wu JY, Weng LH, Li XX, Yu LJ, Xu Y. Valproic acid protects against MPP<sup>+</sup>-mediated neurotoxicity in SH-SY5Y Cells through autophagy. *Neurosci Lett*. 2017 Jan 18;638:60-68
  95. Fan Y, Liu J, Liu D, Zhou Z, Bao Y, Wang J, Zhao Q, Xu Y. NSCA-1-a novel N-substituted coumalamide derivative-increases Adriamycin sensitivity in HepG2/adriamycin cells through modulating Akt/GSK-3  $\beta$  signaling and p53-dependant apoptotic pathway. *ENVIRON TOXICOL CHEM*. 2017 Jan;49:1-7
  96. Chang S, Ruan WC, Xu YZ, Wang YJ, Pang J, Zhang LY, Liao H, Pang T. The natural product 4,10-aromadendranediol induces neuritogenesis in neuronal cells in vitro through activation of the ERK pathway. *Acta Pharmacol Sin*. 2017 Jan;38(1):29-40
  97. Wang Y1, Qiu Y1, Yin S1, Zhang L1, Shi K1, Gao H1, Zhang Z1, He Q1. A functional nanocarrier that copenetrates extracellular matrix and multiple layers of tumor cells for sequential and deep tumor autophagy inhibitor and chemotherapeutic delivery. *Autophagy*. 2017 Feb;13(2):359-370
  98. Wen Y, He W, Jiang M, Zeng M, Cai L. Deriving cells expressing markers of female germ cells from premature ovarian failure patient-specific induced pluripotent stem cells. *Regen Med*. 2017 Mar;12(2):143-152
  99. Zhai Y1, Li Y2, Wang Y3, Cui J2, Feng K4, Kong X2, Chen L5. Psoralidin, a prenylated coumestan, as a novel anti-osteoporosis candidate to enhance boneformation of osteoblasts and decrease bone resorption of osteoclasts. *Eur J Pharmacol*. 2017 Apr 15;801:62-71
  100. Wu J, Yang X, Lu H, Liu L, Xu B, Zheng S, Yu B, Jie K, Wan F. Inhibitory effect and mechanism of exogenous mammalian sterile 20-like kinase 1 on the growth of human colorectal cancer. *Oncol Lett*. 2017 Apr;13(4):2656-2664
  101. Guo Y, Song Z, Zhou M, Yang Y, Zhao Y, Liu B, Zhang X. Infiltrating macrophages in diabetic nephropathy promote podocytes apoptosis via TNF- $\alpha$ -ROS-p38MAPK pathway. *ONCOTARGET*. 2017 Jun 7;8(32):53276-53287
  102. Sun L, Li Y, Jia X, Wang Q, Li Y, Hu M, Tian L, Yang J, Xing W, Zhang W, Wang J, Xu H, Wang L, Zhang D, Ren H. Neuroprotection by IFN- $\gamma$  via astrocyte-secreted IL-6 in acute neuroinflammation. *ONCOTARGET*. 2017 Jun 20;8(25):40065-40078
  103. Xiong LL, Liu F, Lu BT, Zhao WL, Dong XJ, Liu J, Zhang RP, Zhang P, Wang TH. Bone Marrow Mesenchymal Stem-Cell Transplantation Promotes Functional Improvement Associated with CNTF-STAT3 Activation after Hemi-Sectioned Spinal Cord Injury in Tree Shrews. *Front Cell Neurosci*. 2017 Jun 28;11:172.
  104. Wu X, Gao H, Sun W, Yu J, Hu H, Xu Q, Chen X. Nepetoidin B, a Natural Product, Inhibits LPS-stimulated Nitric Oxide Production via Modulation of iNOS Mediated by NF- $\kappa$ B/MKP-5 Pathways. *Phytther Res*. 2017 Jul;31(7):1072-1077
  105. Yan X, Xun M, Dou X, Wu L, Han Y, Zheng J. Regulation of Na<sup>+</sup>-K<sup>+</sup>-ATPase effected high glucose-induced myocardial cell injury through c-Src-dependent NADPH oxidase/ROS pathway. *Exp Cell Res*. 2017 Aug 15;357(2):243-251
  106. Liu JC, Lai FN, Li L, Sun XF, Cheng SF, Ge W, Wang YF, Li L, Zhang XF, De Felici M, Dyce PW, Shen W. Di (2-ethylhexyl) phthalate exposure impairs meiotic progression and DNA damage repair in fetal mouse oocytes in vitro. *Cell Death Dis*. 2017 Aug 3;8(8):e2966
  107. Liu XY, Zhang LJ, Chen Z, Liu LB. The PTEN inhibitor bpV(pic) promotes neuroprotection against amyloid  $\beta$ -peptide(25-35)-induced oxidative stress and neurotoxicity. *Neurol Res*. 2017 Aug;39(8):758-765
  108. Shan ZL, Zhong L, Xiao CL, Gan LG, Xu T, Song H, Yang R, Li L, Liu BZ. Shikonin suppresses proliferation and induces apoptosis in human leukemia NB4 cells through modulation of MAPKs and c-Myc. *Mol Med Rep*. 2017 Sep;16(3):3055-3060
  109. Hao J, Yang Z, Wang L, Zhang Y, Shu Y, Jiang L, Hu Y, Lv W, Dong P, Liu Y. Downregulation of BRD4 inhibits gallbladder cancer proliferation and metastasis and induces apoptosis via PI3K/AKT pathway. *Int J Oncol*. 2017 Sep;51(3):823-831
  110. Liu X, Qiu Z, Wang Z, Zuo W, Gong Z, Liu C, Zeng Q, Qian Y, Jiang L, Li Y, Bu Y, Hu G. NFBD1/MDC1 participates in the regulation of proliferation and apoptosis in human laryngeal squamous cell carcinoma. *Clin Transl Oncol*. 2017 Sep 18
  111. Ying C, Xiao BD, Qin Y, Wang BR, Liu XY, Wang RW, Fang L, Yan H, Zhou XM, Wang YG. GOLPH2-regulated oncolytic adenovirus, GD55, exerts strong killing effect on human prostate cancer stem-like cells in vitro and in vivo. *Acta Pharmacol Sin*. 2017 Sep 7
  112. Liang J, Wang WF, Xie S, Zhang XL, Qi WF, Zhou XP, Hu JX, Shi Q, Yu RT.  $\beta$ -transducin repeat-containing E3 ubiquitin protein ligase inhibits migration, invasion and proliferation of glioma cells. *Oncol Lett*. 2017 Sep;14(3):3131-3135
  113. Geng YD, Zhang C, Lei JL, Yu P, Xia YZ, Zhang H, Yang L, Kong LY. Walsuronoid B induces mitochondrial and lysosomal dysfunction leading to apoptotic rather than autophagic cell death via ROS/p53 signaling pathways in liver cancer. *Biochem Pharmacol*. 2017 Oct 15;142:71-86
  114. Chen GH, Xu CS, Zhang J, Li Q, Cui HH, Li XD, Chang LP, Tang RJ, Xu JY, Tian XQ, Huang PS, Xu J, Jin C, Yang YJ. Inhibition of miR-128-3p by Tongxinluo Protects Human Cardiomyocytes from Ischemia/reperfusion Injury via Upregulation of p70s6k1/p70s6k1. *Front Pharmacol*. 2017 Oct 30;8:775
  115. Liu W, Wang X, Liu Z, Wang Y, Yin B, Yu P, Duan X, Liao Z, Chen Y, Liu C, Li X, Dai Y, Tao Z. SGK1 inhibition induces autophagy-dependent apoptosis via the mTOR-Foxo3a pathway. *BRIT J CANCER*. 2017 Oct 10;117(8):1139-1153
  116. Li Y, Sun Y, Cao T, Su Q, Li Z, Huang M, Ouyang R, Chang H, Zhang S, Miao Y. A cation-exchange controlled core-shell MnS@Bi2S3 theranostic platform for multimodal imaging guided radiation therapy with hyperthermia boost. *Nanoscale*. 2017 Oct 5;9(38):14364-14375
  117. Qian L, Li X, Ye P, Wang G, Dai W, Liu Y, Gao Q, Shen G. Oxymatrine induces apoptosis and inhibits invasion in Gallbladder carcinoma via PTEN/PI3K/AKT pathway. *Cytotechnology*. 2017 Nov 23
  118. Zhao C, Yao X, Chen X, Wu W, Xi F, Yang G, Yu T. Knockdown of

- ubiquitin D inhibits adipogenesis during the differentiation of porcine intramuscular and subcutaneous preadipocytes. *CELL PROLIFERAT*. 2017 Nov 23
- 119.Yang LL1, Zhao Y1, Luo SM1, Ma JY1, Ge ZJ1, Shen W1, Yin S2.Toxic Effects and Possible Mechanisms of HydrogenSulfide and/or Ammonia on Porcine OocyteMaturation in vitro. *Toxicol Lett*. 2017 Dec 29. pii: S0378-4274(17)31529-1
- 120.Xiao B, Qin Y, Ying C, Ma B, Wang B, Long F, Wang R, Fang L, Wang Y.Combination of oncolytic adenovirus and luteolin exerts synergistic antitumor effects in colorectal cancer cells and a mouse model. *Mol Med Rep*. 2017 Dec;16(6):9375-9382
- 121.Li X, Yan X, Wang Y, Wang J, Zhou F, Wang H, Xie W, Kong H.NLRP3 inflammasome inhibition attenuates silica-induced epithelial to mesenchymal transition (EMT) in human bronchial epithelial cells. *Exp Cell Res*. 2017 Dec 16. pii: S0014-4827(17)30668-7
- 122.Xu Z, Zheng M, Zhang L, Zhang X, Zhang Y, Cui X, Gong X, Xi R, Bai R.Effect of ATP and Bax on the apoptosis of *Eimeria tenella* host cells. *BMC Vet Res*. 2017 Dec 28;13(1):399
- 123.Zhang TY,Wu RY,Zhao Y,Xu CS,Zhang WD,Ge W,Liu J,Sun ZY,Zou SH,Shen W.Ochratoxin A exposure decreased sperm motility via the AMPK and PTEN signaling pathways. *TOXICOL APPL PHARM*. 2018 Feb 1;340:49-57
- 124.Sun Y, Wang C, Meng Q, Liu Z, Huo X, Sun P, Sun H, Ma X, Peng J, Liu K..Targeting P-glycoprotein and SORCIN: Dihydromyricetin strengthens anti-proliferative efficiency of adriamycin via MAPK/ERK and Ca<sup>2+</sup>-mediated apoptosis pathways in MCF-7/ADR and K562/ADR. *J Cell Physiol*. 2018 Apr;233(4):3066-3079
- 125.Wang JJ,Yu XW,Wu RY,Sun XF,Cheng SF,Ge W,Liu JC,Li YP,Liu J,Zou SH,De Felici M,Shen W.Starvation during pregnancy impairs fetal oogenesis and folliculogenesis in offspring in the mouse. *Cell Death Dis*. 2018 May 1;9(5):452
- 126.Yi WEI,Xiang-Liang T,Yu Z,Bin L,Lian-Ju S,Chun-Lan L,Tao LIN,Da-Wei HE,Sheng-de WU,Guang-Hui WEI.DEHP exposure destroys blood-testis barrier (BTB) integrity of immature testes through excessive ROS-mediated autophagy. *Genes & Diseases*. 2018 Jun 25;5(3):263-274.
- 127.Tang X,Wu S,Shen L,Wei Y,Cao X,Wang Y,Long C,Zhou Y,Li D,Huang F,Liu B,Wei G.Di-(2-ethylhexyl) phthalate (DEHP)-induced testicular toxicity through Nrf2-mediated Notch1 signaling pathway in Sprague-Dawley rats. *Environ Toxicol*. 2018 Jul;33(7):720-728
- 128.Zhang GL,Song JL,Zhou Y,Zhang RQ,Cheng SF,Sun XF,Qin GQ,Shen W,Li L.Differentiation of sow and mouse ovarian granulosa cells exposed to zearalenone in vitro using RNA-seq gene expression. *TOXICOL APPL PHARM*. 2018 Jul 1;350:78-90
- 129.Zhang GL,Song JL,Ji CL,Feng YL,Yu J,Nyachoti CM,Yang GS.Zearalenone Exposure Enhanced the Expression of Tumorigenesis Genes in Donkey Granulosa Cells via the PTEN/PI3K/AKT Signaling Pathway. *Front Genet*. 2018 Jul 31;9:293
- 130.Zhai QY,Ge W,Wang JJ,Sun XF,Ma JM,Liu JC,Zhao Y,Feng YZ,Dyce PW,De Felici M,Shen W.Exposure to Zinc oxide nanoparticles during pregnancy induces oocyte DNA damage and affects ovarian reserve of mouse offspring. *AGING-US*. 2018 Aug 28;10(8):2170-2189
- 131.Zhang T,Zhou Y,Li L,Zhao Y,De Felici M,Reiter RJ,Shen W.Melatonin protects prepuberal testis from deleterious effects of bisphenol A or diethylhexyl phthalate by preserving H3K9 methylation. *J Pineal Res*. 2018 Sep;65(2):e12497
- 132.Shen L,Tang X,Wei Y,Long C,Tan B,Wu S,Sun M,Zhou Y,Cao X,Wei G.Vitamin E and vitamin C attenuate Di-(2-ethylhexyl) phthalate-induced blood-testis barrier disruption by p38 MAPK in immature SD rats. *Reprod Toxicol*. 2018 Oct;81:17-27
- 133.Zhang RQ,Sun XF,Wu RY,Cheng SF,Zhang GL,Zhai QY,Liu XL,Zhao Y,Shen W,Li L.Zearalenone exposure elevated the expression of tumorigenesis genes in mouse ovarian granulosa cells. *TOXICOL APPL PHARM*. 2018 Oct 1;356:191-203
- 134.Zhao Y,Zhao R,Wu J,Wang Q,Pang K,Shi Q,Gao Q,Hu Y,Dong X,Zhang J,Sun J.Melatonin protects against A $\beta$  -induced neurotoxicity in primary neurons via miR-132/PTEN/AKT/FOXO3a pathway. *Biofactors*. 2018 Nov;44(6):609-618
- 135.Zhou B,Zhang JY,Liu XS,Chen HZ,Ai YL,Cheng K,Sun RY,Zhou D,Han J,Wu Q.Tom20 senses iron-activated ROS signaling to promote melanoma cell pyroptosis. *Cell Res*. 2018 Dec;28(12):1171-1185
- 136.Sun ZY,Zhang P,Wang JJ,Liu JC,Li L,Shen W,Zhai QY.Melatonin alleviates meiotic defects in fetal mouse oocytes induced by Di (2-ethylhexyl) phthalate in vitro. *AGING-US*. 2018 Dec 26;10(12):4175-4185
- 4187
- 137.Zhu P,Liu Z,Zhou J,Chen Y.Tanshinol inhibits the growth, migration and invasion of hepatocellular carcinoma cells via regulating the PI3K-AKT signaling pathway. *ONCOTARGETS THER*. 2018 Dec 19;12:87-99
- 138.Yang LL,Cui YX,Ma JY,Ge ZJ,Shen W,Yin S.Tributyltin oxide exposure impairs mouse oocyte maturation and its possible mechanisms. *J Cell Biochem*. 2019 Jan;120(1):715-726
- 139.Chu P,He L,Zhu D,Chen L,Huang R,Liao L,Li Y,Zhu Z,Wang Y.Identification, characterisation and preliminary functional analysis of IRAK-M in grass carp (*Ctenopharyngodon idella*). *FISH SHELLFISH IMMUN*. 2019 Jan;84:312-321
- 140.Zhou R,Wang H,Yang Y,Zhang C,Dong X,Du J,Yan L,Zhang G,Gu Z,Zhao Y.Tumor microenvironment-manipulated radiocatalytic sensitizer based on bismuth heteropolytungstate for radiotherapy enhancement. *Biomaterials*. 2019 Jan;189:11-22
- 141.Han B,Xu K,Liu Z,Ge W,Shao S,Li P,Yan N,Li X,Zhang Z.Oral yeast-based DNA vaccine confers effective protection from *Aeromonas hydrophila* infection on *Carassius auratus*. *FISH SHELLFISH IMMUN*. 2019 Jan;84:948-954
- 142.Su W,Wang Y,Wang F,Zhang B,Zhang H,Shen Y,Yang H.Circular RNA hsa\_circ\_0007059 indicates prognosis and influences malignant behavior via AKT/mTOR in oral squamous cell carcinoma. *J Cell Physiol*. 2019 Jan 24
- 143.Yang R,Xing L,Zheng X,Sun Y,Wang X,Chen J.The circRNA circAGFG1 acts as a sponge of miR-195-5p to promote triple-negative breast cancer progression through regulating CCNE1 expression. *Mol Cancer*. 2019 Jan 8;18(1):4
- 144.Jiang Y,Zhang T,Vallesi A,Yang X,Gao F.Time-course analysis of nuclear events during conjugation in the marine ciliate *Euploites vannus* and comparison with other ciliates (Protozoa, Ciliophora). *Cell Cycle*. 2019 Feb;18(3):288-298
- 145.Cao P,Feng Y,Deng M,Li J,Cai H,Meng Q,Fang W,Li Y,Ke J,Long X.MiR-15b is a key regulator of proliferation and apoptosis of chondrocytes from patients with condylar hyperplasia by targeting IGF1, IGF1R and BCL2. *OSTEOARTHR CARTILAGE*. 2019 Feb;27(2):336-346
- 146.Liang L,Hui K,Hu C,Wen Y,Yang S,Zhu P,Wang L,Xia Y,Qiao Y,Sun W,Fei J,Chen T,Zhao F,Yang B,Jiang X.Autophagy inhibition potentiates the anti-angiogenic property of multi-kinase inhibitor anlotinib through JAK2/STAT3/VEGFA signaling in non-small cell lung cancer cells. *J EXP CLIN CANC RES*. 2019 Feb 12;38(1):71
- 147.Zheng Y,Tu C,Zhang J,Wang J.Inhibition of multiple myeloma-derived exosomes uptake suppresses the functional response in bone marrow stromal cell. *Int J Oncol*. 2019 Mar;54(3):1061-1070
- 148.Ruan W,Li J,Xu Y,Wang Y,Zhao F,Yang X,Jiang H,Zhang L,Saavedra JM,Shi L,Pang T.MALAT1 Up-Regulator Polydatin Protects Brain Microvascular Integrity and Ameliorates Stroke Through C/EBP $\beta$ /MALAT1/CREB/PGC-1 $\alpha$ /PPAR $\gamma$  Pathway. *Cell Mol Neurobiol*. 2019 Mar;39(2):265-286
- 149.Tian J,Rui K,Hong Y,Wang X,Xiao F,Lin X,Ma J,Guo H,Xu H,Ma K,Xu D,Liu D,Zhao Y,Lu L,Wang S.Increased GITRL Impairs the Function of Myeloid-Derived Suppressor Cells and Exacerbates Primary Sjögren Syndrome. *J Immunol*. 2019 Mar 15;202(6):1693-1703
- 150.Sun Y,Liu W,Wang C,Meng Q,Liu Z,Huo X,Yang X,Sun P,Sun H,Ma X,Peng J,Liu K.Combination of dihydromyricetin and ondansetron strengthens antiproliferative efficiency of adriamycin in K562/ADR through downregulation of SORCIN: A new strategy of inhibiting P-glycoprotein. *J Cell Physiol*. 2019 Apr;234(4):3685-3696
- 151.Huang Z,Li Q,Yuan Y,Zhang C,Wu L,Liu X,Cao W,Guo H,Duan S,Xu X,Zhang B,Xing C.Renalase attenuates mitochondrial fission in cisplatin-induced acute kidney injury via modulating sirtuin-3. *Life Sci*. 2019 Apr 1;222:78-87
- 152.Su W,Wang Y,Wang F,Sun S,Li M,Shen Y,Yang H.Hsa\_circ\_0005379 regulates malignant behavior of oral squamous cell carcinoma through the EGFR pathway. *BMC Cancer*. 2019 Apr 29;19(1):400
- 153.Zhu D,Huang R,Fu P,Chen L,Luo L,Chu P,He L,Li Y,Liao L,Zhu Z,Wang Y.Investigating the Role of BATF3 in Grass Carp (*Ctenopharyngodon idella*) Immune Modulation: A Fundamental Functional Analysis. *Int J Mol Sci*. 2019 Apr 4;20(7). pii: E1687
- 154.Li WD,Yu S,Luo SM,Shen W,Yin S,Sun QY.Melatonin defends mouse oocyte quality from benzo[ghi]perylene-induced deterioration. *J Cell Physiol*. 2019 May;234(5):6220-6229
- 155.Chen C,Liu Q,Hua H,Wang X,Wang P,Cui Z,Qian T.Novel microRNA,

- miR-sc6, modulates Schwann cell phenotype via targeting ErbB4.*Exp Ther Med.* 2019 May;17(5):4116-4122
- 156.Zhou X,Li T,Chen Y,Zhang N,Wang P,Liang Y,Long M,Liu H,Mao J,Liu Q,Sun X,Chen H.Mesenchymal stem cell-derived extracellular vesicles promote the in vitro proliferation and migration of breast cancer cells through the activation of the ERK pathway.*Int J Oncol.* 2019 May;54(5):1843-1852
- 157.Chen C,Ai Q,Chu S,Zhang Z,Zhou X,Luo P,Liu Y,Chen N.IMM-H004 protects against oxygen-glucose deprivation/reperfusion injury to BV2 microglia partly by modulating CKLF1 involved in microglia polarization.*Int Immunopharmacol.* 2019 May;70:69-79
- 158.Li J,Xing F,Chen F,He L,So KF,Liu Y,Xiao J.Functional 3D Human Liver Bud Assembled from MSC-Derived Multiple Liver Cell Lineages.*Cell Transplant.* 2019 May;28(5):510-521
- 159.Liu Y,Bi T,Liu L,Gao Q,Shen G,Qin L.S-Adenosylmethionine synergistically enhances the antitumor effect of gemcitabine against pancreatic cancer through JAK2/STAT3 pathway.*N-S ARCH PHARMACOL.* 2019 May;392(5):615-622
- 160.Chu P,He L,Yang C,Li Y,Huang R,Liao L,Li Y,Zhu Z,Wang Y.Characterisation and function of TRIM23 in grass carp (*Ctenopharyngodon idella*).*FISH SHELLFISH IMMUN.* 2019 May;88:627-635
- 161.Liu H,Zhao J,Fu R,Zhu C,Fan D.The ginsenoside Rk3 exerts anti-esophageal cancer activity in vitro and in vivo by mediating apoptosis and autophagy through regulation of the PI3K/Akt/mTOR pathway.*PLoS One.* 2019 May 15;14(5):e0216759
- 162.Ye M,Zhu X,Yan R,Hui J,Zhang J,Liu Y,Li X.Sodium demethylcantharidate induces apoptosis in hepatocellular carcinoma cells via ER stress.*Am J Transl Res.* 2019 May 15;11(5):3150-3158.eCollection 2019
- 163.Qimuge N,He Z,Qin J,Sun Y,Wang X,Yu T,Dong W,Yang G,Pang W.Overexpression of DNMT3A promotes proliferation and inhibits differentiation of porcine intramuscular preadipocytes by methylating p21 and PPAR $\gamma$  promoters.*Gene.* 2019 May 15;696:54-62
- 164.Yan H,Ma L,Wang H,Wu S,Huang H,Gu Z,Jiang J,Li Y.Luteolin decreases the yield of influenza A virus in vitro by interfering with the coat protein I complex expression.*J NAT MED-TOKYO.* 2019 Jun;73(3):487-496
- 165.Shen Z,Chen J,Lu H.Osthole induced apoptosis in human normal liver cells by regulating cell proliferation and endoplasmic reticulum stress.*Environ Toxicol.* 2019 Jun;34(6):768-776
- 166.Zhang S,Jiang H,Huang S,Li P,Wang F,Curdian sulfate/O-linked quaternary chitosan nanoparticles acting as potential adjuvants promote multiple arms of immune responses.*CARBOHYD POLYM.* 2019 Jun 1;213:100-111
- 167.Tang L,Lu W,Huang J,Tang X,Zhang H,Liu S.miR-144 promotes the proliferation and differentiation of bone mesenchymal stem cells by downregulating the expression of SFRP1.*Mol Med Rep.* 2019 Jul;20(1):270-280
- 168.Liu S,Xing Y,Wang J,Pan R,Li G,Tang H,Chen G,Yan L,Guo L,Jiang M,Gong Z,Lin L,Dong J.The Dual Role of HIV-1 gp120 V3 Loop-Induced Autophagy in the Survival and Apoptosis of the Primary Rat Hippocampal Neurons.*Neurochem Res.* 2019 Jul;44(7):1636-1652
- 169.Zhu D,Li Y,Huang R,Luo L,Chen L,Fu P,He L,Li Y,Liao L,Zhu Z,Wang Y,Molecular characterization and functional activity of Prx1 in grass carp (*Ctenopharyngodon idella*).*FISH SHELLFISH IMMUN.* 2019 Jul;90:395-403
- 170.Zhu J,Ye Q,Xu S,Chang YX,Liu X,Ma Y,Zhu Y,Hua S.Shengmai injection alleviates H2O2-induced oxidative stress through activation of AKT and inhibition of ERK pathways in neonatal rat cardiomyocytes.*J Ethnopharmacol.* 2019 Jul 15;239:111677
- 171.Wu S,Wang HQ,Guo TT,Li YH.Luteolin inhibits CVB3 replication through inhibiting inflammation.*J Asian Nat Prod Res.* 2019 Jul 19:1-12
- 172.Tan T,Hu H,Wang H,Li J,Wang Z,Wang J,Wang S,Zhang Z,Li Y.Bioinspired lipoproteins-mediated photothermia remodels tumor stroma to improve cancer cell accessibility of second nanoparticles.*Nat Commun.* 2019 Jul 25;10(1):3322
- 173.Zhang CY,Gao Y,Zhu RX,Qiao YN,Zhou JC,Zhang JZ,Li Y,Li SW,Fan SH,Lou HX.Prenylated Bibenzyls from the Chinese Liverwort Radula constricta and Their Mitochondria-Derived Paraptotic Cytotoxic Activities.*J Nat Prod.* 2019 Jul 26;82(7):1741-1751
- 174.Deng JL,Zhang R,Zeng Y,Zhu YS,Wang G.Statins induce cell apoptosis through a modulation of AKT/FOXO1 pathway in prostate cancer cells.*Cancer Manag Res.* 2019 Jul 31;11:7231-7242
- 175.Wang JJ,Zhai QY,Zhang RQ,Ge W,Liu JC,Li L,Sun ZY,De Felici M,Shen W.Effects of activin A on the transcriptome of mouse oogenesis in vitro.*J Cell Physiol.* 2019 Aug;234(8):14339-14350
- 176.Yin W,Yang L,Kong D,Nie Y,Liang Y,Teng CB.Guanine-rich RNA binding protein GRSF1 inhibits myoblast differentiation through repressing mitochondrial ROS production.*Exp Cell Res.* 2019 Aug 1;381(1):139-149
- 177.Chen J,Chen X,Yao J,Li M,Yang X.The combination of Decitabine and EPZ-6438 effectively facilitate adipogenic differentiation of induced pluripotent stem cell-derived mesenchymal stem cells.*BIOCHEM BIOPH RES CO.* 2019 Aug 13;516(1):307-312
- 178.Zhang K,Li H,Dong S,Liu Y,Wang D,Liu H,Su F,Ge L,Jiang Y.Establishment and evaluation of a PRRSV-sensitive porcine endometrial epithelial cell line by transfecting SV40 large T antigen.*BMC Vet Res.* 2019 Aug 19;15(1):299
- 179.Jiang S,Lyu C,Zhao P,Li W,Kong W,Huang C,Genin GM,Du Y.Cryoprotectant enables structural control of porous scaffolds for exploration of cellular mechano-responsiveness in 3D.*Nat Commun.* 2019 Aug 2;10(1):3491
- 180.Duan H,Lei Z,Xu F,Pan T,Lu D,Ding P,Zhu C,Pan C,Zhang S.PARK2 Suppresses Proliferation and Tumorigenicity in Non-small Cell Lung Cancer.*Front Oncol.* 2019 Aug 23;9:790
- 181.Yan-Yun Pan,Jin-Xiu Yang,Wei Mao,Xing-Xiang Wang.RNA-binding protein SFPQ cooperates with HDAC1 to suppress CD40 transcription in pulmonary adventitial fibroblasts.*Cell Biol Int.* 2019 Aug 8.;doi: 10.1002/cbin.11216
- 182.Chu P,He L,Yang C,Zeng W,Huang R,Liao L,Li Y,Zhu Z,Wang Y.Grass carp ATG5 and ATG12 promote autophagy but down-regulate the transcriptional expression levels of IFN- $\gamma$  signaling pathway.*FISH SHELLFISH IMMUN.* 2019 Sep;92:600-611
- 183.Zhang T,He L,Sun W,Qin Y,Zhang P,Zhang H,1,25-Dihydroxyvitamin D3 enhances the susceptibility of anaplastic thyroid cancer cells to adriamycin-induced apoptosis by increasing the generation of reactive oxygen species.*Mol Med Rep.* 2019 Sep;20(3):2641-2648
- 184.Chen R,Hou R,Hong X,Yan S,Zha J.Organophosphate flame retardants (OPFRs) induce genotoxicity in vivo: A survey on apoptosis, DNA methylation, DNA oxidative damage, liver metabolites, and transcriptomics.*Environ Int.* 2019 Sep;130:104914
- 185.Hou DD,Zhang W,Gao YL,Sun YZ,Wang HX,Qi RQ,Chen HD,Gao XH.Anti-inflammatory effects of quercetin in a mouse model of MC903-induced atopic dermatitis.*Immunopharmacol.* 2019 Sep;74:105676
- 186.Zhu D,Huang R,Yang C,Fu P,Chen L,Jiang Y,He L,Li Y,Liao L,Zhu Z,Wang Y.Identification and molecular characterization of peroxiredoxin 2 in grass carp (*Ctenopharyngodon idella*).*FISH SHELLFISH IMMUN.* 2019 Sep;92:570-582
- 187.Pan JA,Tang Y,Yu JY,Zhang H,Zhang JF,Wang CQ,Gu J.miR-146a attenuates apoptosis and modulates autophagy by targeting TAF9b/P53 pathway in doxorubicin-induced cardiotoxicity.*Cell Death Dis.* 2019 Sep 11;10(9):668
- 188.Zhong B,Shi D,Wu F,Wang S,Hu H,Cheng C,Qing X,Huang X,Luo X,Zhang Z,Shao Z.Dynasore suppresses cell proliferation, migration, and invasion and enhances the antitumor capacity of cisplatin via STAT3 pathway in osteosarcoma.*Cell Death Dis.* 2019 Sep 18;10(10):687
- 189.Li Y,Ruan X,Xu X,Li C,Qiang T,Zhou H,Gao J,Wang X.Shengmai Injection Suppresses Angiotensin II-Induced Cardiomyocyte Hypertrophy and Apoptosis via Activation of the AMPK Signaling Pathway Through Energy-Dependent Mechanisms.*Front Pharmacol.* 2019 Sep 20;10:1095
- 190.Zhan W,Liao X,Li L,Chen Z,Tian T,Yu L,Chen Z.In vitro mitochondrial-targeted antioxidant peptide induces apoptosis in cancer cells.*ONCOTARGETS THER.* 2019 Sep 6;12:7297-7306
- 191.Yang L,Zhu Y,Kong D,Gong J,Yu W,Liang Y,Nie Y,Teng CB.EGF suppresses the expression of miR-124a in pancreatic  $\beta$  cell lines via ETS2 activation through the MEK and PI3K signaling pathways.*Int J Biol Sci.* 2019 Sep 7;15(12):2561-2575
- 192.Chen R,Huang H,Liu H,Xi J,Ning J,Zeng W,Shen C,Zhang T,Yu G,Xu Q,Chen X,Wang J,Lu F.Friend or Foe? Evidence Indicates Endogenous Exosomes Can Deliver Functional gRNA and Cas9 Protein.*Small.* 2019 Sep;15(38):e1902686
- 193.Jing ZF,Bi JB,Li Z,Liu X,Li J,Zhu Y,Zhang XT,Zhang Z,Li Z,Kong CZ.Inhibition of miR-34a-5p can rescue disruption of the p53-DAPK axis to suppress progression of clear cell renal cell carcinoma.*Mol*

194. Wang D,Du Y,Xu H,Pan H,Wang R.Paeonol protects mitochondrial injury and prevents pulmonary vascular remodeling in hypoxia.RESP PHYSIOL NEUROBI. 2019 Oct;268:103252
195. Li J,Xu S,Liu Y,Yan Z,Zhang F,Lv Q,Tong N.Activated PPAR  $\beta$  /  $\delta$  Protects Pancreatic  $\beta$  Cells in Type 2 Diabetic Goto-Kakizaki Rats from Lipoapoptosis via GPR40.Lipids. 2019 Oct;54(10):603-616
196. Zhang L,Zheng MX,Xi R,Xu ZY,Zhang XS,Zheng LL,Bai R,Mi CL,Hao FF,Feng YP.Comparison of the host cells apoptosis induced by precocious strains and virulent strains of *Eimeria tenella*.J POULT SCI. 2019 Oct 1;98(10):4384-4390
197. Yin L,Chen X,Ji H,Gao S.Dexmedetomidine protects against sepsis-associated encephalopathy through Hsp90/AKT signaling.Mol Med Rep. 2019 Nov;20(5):4731-4740
198. Li N,Wang LJ,Xu WL,Liu S,Yu JY.MicroRNA-379-5p suppresses renal fibrosis by regulating the LIN28/let-7 axis in diabetic nephropathy.Int J Mol Med. 2019 Nov;44(5):1619-1628
199. Deng H,Zhang W,Li J,Li J,Hu L,Yan W,Liu S,He J,Weng S.A signal transducers and activators of transcription (STAT) gene from *Scylla paramamosain* is involved in resistance against mud crab reovirus.FISH SHELLFISH IMMUN. 2019 Nov;94:580-591
200. Zhu D,Fu P,Huang R,Xiong L,Wang Y,He L,Liao L,Li Y,Zhu Z,Wang Y.Molecular characterization, tissue distribution and functional analysis of galectin 1-like 2 in grass carp (Ctenopharyngodon idella).FISH SHELLFISH IMMUN. 2019 Nov;94:455-463
201. Fan J,Wang Q,Zhang Z,Sun L.Curcumin mitigates the epithelial-to-mesenchymal transition in biliary epithelial cells through upregulating CD109 expression.Drug Dev Res. 2019 Nov;80(7):992-999.
202. Ying J,Han Z,Pei S,Su L,Ruan D.Effects of Stromal Cell-Derived Factor-1 $\alpha$  Secreted in Degenerative Intervertebral Disc on Activation and Recruitment of Nucleus Pulposus-Derived Stem Cells.Stem Cells Int. 2019 Nov 19;2019:9147835
203. Tang Q, Chen Z, Zhao L, Xu H.Circular RNA hsa\_circ\_0000515 acts as a miR-326 sponge to promote cervical cancer progression through up-regulation of ELK1.AGING-US. 2019 Nov 26;11(22):9982-9999
204. Chu P,He L,Zhu D,Huang R,Liao L,Li Y,Zhu Z,Wang Y.Identification, expression and functional characterisation of CYP1A in grass carp (Ctenopharyngodon idella).FISH SHELLFISH IMMUN. 2019 Dec;95:35-43
205. Han X,Wang R,Xu J,Chen Q,Liang C,Chen J,Zhao J,Chu J,Fan Q,Archibong E,Jiang L,Wang C,Liu Z.In situ thermal ablation of tumors in combination with nano-adjuvant and immune checkpoint blockade to inhibit cancer metastasis and recurrence.Biomaterials. 2019 Dec;224:119490
206. Yu X,Hao M,Liu Y,Ma X,Lin W,Xu Q,Zhou H,Shao N,Kuang H.Liraglutide ameliorates non-alcoholic steatohepatitis by inhibiting NLRP3 inflammasome and pyroptosis activation via mitophagy.Eur J Pharmacol. 2019 Dec 1;864:172715
207. Guo Q,Li J,Zhong L,Cai H,Wang H,Yan H,Xia G,Kang Y,Wang C.Prostaglandin-E2 deficiency during late pregnancy and the associated increase in interleukin-1  $\beta$  derived from periaortic lymph nodes lead to abortion.Mol Hum Reprod. 2019 Dec 1;25(12):825-837
208. Li T,Chen Q,Zheng Y,Zhang P,Chen X,Lu J,Lv Y,Sun S,Zeng W.PAMAM-cRGD mediating efficient siRNA delivery to spermatogonial stem cells.Stem Cell Res Ther. 2019 Dec 18;10(1):399
209. Zhang N,Zeng X,Sun C,Guo H,Wang T,Wei L,Zhang Y,Zhao J,Ma X.LncRNA LINC00963 Promotes Tumorigenesis and Radioresistance in Breast Cancer by Sponging miR-324-3p and Inducing ACK1 Expression.MOL THER-NUCL ACIDS. 2019 Dec 6;18:871-881
210. Li J,Yang X,Liu F,Song Y,Liu Y.Activated PI3K/AKT Signaling Pathway Associates with Oxidative Stress and Impaired Developmental Potential of Vitrified-Thawed Oocytes.Reprod Sci. 2020 Jan;27(1):404-410
211. Xu W,Zhao Y,Pan B,Liu Y,Li Y,Bourland WA,Luo X.Morphology, Morphogenesis, and Phylogeny of Urosoma caudata (Ehrenberg, 1833) Berger, 1999 (Ciliophora, Hypotrichia) based on a Chinese Population.J Eukaryot Microbiol. 2020 Jan;67(1):76-85
212. Zhang B,Zhu L,Dai Y,Li H,Huang K,Luo Y,Xu W.An in vitro attempt at precision toxicology reveals the involvement of DNA methylation alteration in ochratoxin A-induced G0/G1 phase arrest.Epigenetics. 2020 Jan - Feb;15(1-2):199-214
213. Tang C,Ou J,Kou L,Deng J,Luo S.Circ\_016719 plays a critical role in neuron cell apoptosis induced by I/R via targeting miR-29c/Map2k6.MOL CELL PROBE. 2020 Feb;49:101478
214. Zhang Y,Huang WJ,Yang QR,Zhang HD,Zhu XJ,Zeng M,Zhou X,Wang ZY,Li WJ,Jing HS,Zhang XB,Shi YP,Hu H,Yan HX,Li ZH,Zhai B.Cryopreserved biopsy tissues of rectal cancer liver metastasis for assessment of anticancer drug response in vitro and in vivo.Oncol Rep. 2020 Feb;43(2):405-414
215. Yi Li,Jing Lu,Xin Cao,Hongwei Zhao,Longfei Gao,Peng Xia,Gang Pei.A Newly Synthesized Rhamnoside Derivative Alleviates Alzheimer's Amyloid-  $\beta$ -Induced Oxidative Stress, Mitochondrial Dysfunction, and Cell Senescence through Upregulating SIRT3Oxid Med Cell Longev. 2020 Feb 13;2020:7698560.;doi: 10.1155/2020/7698560
216. Zhu D,Huang R,Chu P,Chen L,Li Y,He L,Li Y,Liao L,Zhu Z,Wang Y.Characterization and expression of galectin-3 in grass carp (Ctenopharyngodon idella).Dev Comp Immunol. 2020 Mar;104:103567
217. Wang XL,Ma YX,Xu RJ,Ma JJ,Zhang HC,Qi SB,Xu JH,Qin XZ,Zhang HN,Liu CM,Chen JQ,Li B,Yang HL,Sajilafu.c-Myc controls the fate of neural progenitor cells during cerebral cortex development.J Cell Physiol. 2020 Apr;235(4):4011-4021
218. Siyu Li,Xiaomeng Ren,Yuandong Wang,Jiangning Hu,Haitao Wu,Shuang Song,Chunhong Yan.Fucoxanthin alleviates palmitate-induced inflammation in RAW 264.7 cells through improving lipid metabolism and attenuating mitochondrial dysfunction Food Funct. 2020 Apr 30;11(4):3361-3370.;doi: 10.1039/d0fo00442a
219. Deng H,Hu L,Li J,Yan W,Song E,Kuang M,Liu S,He J,Weng S.The NF-  $\kappa$  B family member dorsal plays a role in immune response against Gram-positive bacterial infection in mud crab (*Scylla paramamosain*).Dev Comp Immunol. 2020 May;106:103581
220. Tian-Yu Zhang,Li Kong,Jia-Xing Hao,Han Wang,Zi-Hui Yan,Xiao-Feng Sun,Wei Shen.Effects of Ochratoxin A exposure on DNA damage in porcine granulosa cells in vitro Toxicol Lett. 2020 May 23;330:167-175.;doi: 10.1016/j.toxlet.2020.05.011
221. Yonglu Li,Yong Jiang,Qiang Chu,Xiaodong Zheng.Radix Tetragastris extract from different origins protect RAW264.7 macrophages against LPS-induced inflammation J Food Sci. 2020 May;85(5):1586-1595.;doi: 10.1111/1750-3841.15113
222. Chunming Zhao,Linlin Li,Zhiwei Li,Jiawen Xu,Qing Yang,Peng Shi,Kaining Zhang,Rui Jiang.A Novel Circular RNA hsa\_circRPPH1\_015 Exerts an Oncogenic Role in Breast Cancer by Impairing miRNA-326-Mediated ELK1 Inhibition Front Oncol. 2020 Jun 24;10:906.;doi: 10.3389/fonc.2020.00906
223. He Ding,Yang Song,Xiaowan Huang,Liansheng Wang,Shanzi Luo,Hao Zhang,Hao Pan,Wenwei Jiang,Jing Qian,Guangyu Yao,Longping Wen,Yunjiao Zhang.mTORC1-dependent TFEB nucleus translocation and pro-survival autophagy induced by zeolitic imidazolate framework-8 BIOMATER SCI-UK. 2020 Jul 28;8(15):4358-4369.;doi: 10.1039/d0bm00773k
224. Yuqing Zhang,Huanfeng Bian,Yu Ma,Yuanyuan Xiao,Fang Xiao,Cr(VI)-induced overactive mitophagy contributes to mitochondrial loss and cytotoxicity in L02 hepatocytes Biochem J. 2020 Jul 31;477(14):2607-2619.;doi: 10.1042/BCJ20200262. (IF 4.097)
225. Wei-Dong Li,Chuan-Jie Zang,Shen Yin,Wei Shen,Qing-Yuan Sun,Minghui Zhao.Metformin protects against mouse oocyte apoptosis defects induced by arecoline CELL PROLIFERAT. 2020 Jul;53(7):e12809.;doi: 10.1111/cpr.12809. (IF 5.753)
226. Xiaochen Qiu,Kaisheng Dong,Jingzhi Guan,JianMiao He.Hydrogen attenuates radiation-induced intestinal damage by reducing oxidative stress and inflammatory response Int Immunopharmacol. 2020 Jul;84:106517.;doi: 10.1016/j.intimp.2020.106517
227. Wen-Xiang Liu,Shao-Jing Tan,Yu-Feng Wang,Lan Li,Xiao-Feng Sun,Jing Liu,Francesca Gioia Klinger,Massimo De Felici,Wei Shen,Shun-Feng Cheng.Melatonin ameliorates murine fetal oocyte meiotic dysfunction in F1 and F2 offspring caused by nicotine exposure during pregnancy Environ Pollut. 2020 Aug;263(Pt A):114519.;doi: 10.1016/j.envpol.2020.114519
228. Qinlin Shi,Huan Wu,Yonglin Li,Lianju Shen,Xiaomao Tian,Tao Lin,Guanghui Wei.Inhibition of Wilms' Tumor Proliferation and Invasion by Blocking TGF-  $\beta$  Receptor I in the TGF-  $\beta$ /Smad Signaling Pathway Biomed Res Int. 2020 Nov 16;2020:8039840.;doi: 10.1155/2020/8039840
229. Jing Huang,Ting Ting Zhang,Ke Jiang,Wan Shu Hong,Shi Xi Chen GFP expression pattern in pituitary and gonads under the control of nuclear progesterone receptor promoter in transgenic zebrafish Dev Dyn. 2020 Nov;249(11):1365-1376.;doi: 10.1002/dvdy.213.
230. Chao Tian,Shuai Li,Lang He,Xiaobo Han,Feng Tang,Rongqi

- Huang,Zuoxian Lin,Sihao Deng,Junjie Xu,Hualin Huang,Huifang Zhao,Zhiyuan Li.Transient receptor potential ankyrin 1 contributes to the lysophosphatidylcholine-induced oxidative stress and cytotoxicity in OLN-93 oligodendrocyte CELL STRESS CHAPERON. 2020 Nov;25(6):955-968.;doi: 10.1007/s12192-020-01131-y
- 231.Jing-Cai Liu,Zi-Hui Yan,Bo Li,Hong-Chen Yan,Massimo De Felici,Wei Shen.Di (2-ethylhexyl) phthalate impairs primordial follicle assembly by increasing PDE3A expression in oocytes Environ Pollut. 2021 Feb 1;270:116088.;doi: 10.1016/j.envpol.2020.116088
- 232.Yuhao Wu,Junke Wang,Tianxin Zhao,Yuexin Wei,Lindong Han,Lianju Shen,Chunlan Long,Shengde Wu,Guanghui Wei.LncRNAs activate longevity regulation pathway due to aging of Leydig cells caused by DEHP exposure: A transcriptome-based study ECOTOX ENVIRON SAFE. 2021 Feb;209:111798.;doi: 10.1016/j.ecoenv.2020.111798
- 233.Wenjun Yu,Busong Wang,Liang Zhou,Guoqiang Xu.Endoplasmic Reticulum Stress-Mediated p62 Downregulation Inhibits Apoptosis via c-Jun Upregulation BIOMOL THER. 2021 Mar 1;29(2):195-204.;doi: 10.4062/biomolther.2020.089
- 234.Tingting Song,Weiren Yang,Libo Huang,Zaibin Yang,Shuzhen Jiang.Zearalenone exposure affects the Wnt/ $\beta$ -catenin signaling pathway and related genes of porcine endometrial epithelial cells in vitro Anim Biosci. 2021 Jun;34(6):993-1005.;doi: 10.5713/ajas.20.0292.

Version 2021.09.01