碧云天生物技术/Beyotime Biotechnology  
 订货热线: 400-1683301 或 800-8283301  
 订货 e-mail: order@beyotime.com  
 技术咨询: info@beyotime.com  
 网址: http://www.beyotime.com

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1. Wang F, Liang YJ, Wu XP, Su XD, Fu LW. . ABCG2-overexpressing S1-M1-80 cell xenografts in nude mice keep original biochemistry and cellbiological properties. *Chin J Cancer*. 2012 Mar;31(3):150-8.
2. Sun J, Liu X, Gao H, Zhang L, Ji Q, Wang Z, Zhou L, Wang Y, Sui H, Fan Z, Li Q. . Overexpression of colorectal cancer oncogene CHRDL2 predicts a poor prognosis. *ONCOTARGET*. 2017 Feb 14;8(7):11489-11506.
3. Yu DJ, Wang XJ, Shi YF, Jiang CY, Zhao RZ, Zhu YP, Chen L, Yang YQ, Sun XW, Xia SJ. . Macrophages are targets of retinoic acid signaling during the wound-healing process afterthulium laser resection of the prostate. *ONCOTARGET*. 2017 May 26;8(42):71996-72007.
4. Xudong Wang, Taiqiu Chen, Zhihui Deng, Wenjie Gao, Tongzhou Liang, Xianjian Qiu, Bo Gao, Zizhao Wu, Jincheng Qiu, Yuanxin Zhu, Yanbo Chen, Zhancheng Liang, Hang Zhou, Caixia Xu, Anjing Liang, Peiqiang Su, Yan Peng, Dongsheng Huang . Melatonin promotes bone marrow mesenchymal stem cell osteogenic differentiation and prevents osteoporosis development through modulating circ\_0003865 that sponges miR-3653-3p *Stem Cell Res Ther*. 2021 Feb 25;12(1):150.
5. Wenwen Cui, Yuanyuan Hao, Mingye Wang, Qiuyan Zhang, Junmei Wang, Gang Wei, Yunlong Hou . Inhibition of Autophagy Facilitates XY03-EA-Mediated Neuroprotection against the Cerebral Ischemia/Reperfusion Injury in Rats *Oxid Med Cell Longev*. 2022 Mar 30;2022:7013299.
6. Nie J, Cao X, Zhou M, Zhang X, Zhang R, Niu L, Xia Y, Hou L, Wang C. . Musca domestica pupae lectin induces apoptosis in HepG2 cells through a NF- $\kappa$ B/p65-caspase pathway. *B CANCER*. 2012 Apr 1;99(4):E49-54.
7. Meng X, Li F, Chen S, Tang C, Zhang W, Wang Z, Zhao S. . Cloning and expression of neuron-specific enolase in the corpus luteum of dairy goats. *Gene*. 2012 Jul 25;503(2):222-8.
8. Zhang W1, Wu Q1, Lu YL1, Gong QH1, Zhang F1, Shi JS1. . Protective effects of *Dendrobium nobile* Lindl. alkaloids on amyloid beta (25-35)-induced neuronal injury. *Neural Regen Res*. 2017 Jul;12(7):1131-1136.
9. Fan DX, Yang XH, Li YN, Guo L . 17 $\beta$ -Estradiol on the Expression of G-Protein Coupled Estrogen Receptor (GPER/GPR30) Mitophagy, and the PI3K/Akt Signaling Pathway in ATDC5 Chondrocytes In Vitro. *MED SCI MONITOR*. 2018 Apr 2;24:1936-1947.
10. Dong-Xiao Fan, Xu-Hao Yang, Yi-Nan Li, Lei Guo . 17 $\beta$ -Estradiol on the Expression of G-Protein Coupled Estrogen Receptor (GPER/GPR30) Mitophagy, and the PI3K/Akt Signaling Pathway in ATDC5 Chondrocytes In Vitro *MED SCI MONITOR*. 2018 Apr 2;24:1936-1947.
11. Zhang L, Zhao Y, Guo L. . 17 $\beta$ -estradiol protects INS-1 insulinoma cells from mitophagy via G protein-coupled estrogen receptors and the PI3K/Akt signaling pathway. *Int J Mol Med*. 2018 May;41(5):2839-2846.
12. Zhang L, Zhao Y, Guo L . 17 $\beta$ -estradiol protects INS-1 insulinoma cells from mitophagy via G protein-coupled estrogen receptors and the PI3K/Akt signaling pathway. *Int J Mol Med*. 2018 May;41(5):2839-2846.
13. Shen J, Tu L, Chen D, Tan T, Wang Y, Wang S . TRPV4 channels stimulate Ca<sup>2+</sup>-induced Ca<sup>2+</sup> release in mouse neurons and trigger endoplasmic reticulum stress after intracerebral hemorrhage. *Brain Res Bull*. 2019 Mar 146:143-152.
14. Zhong X, Xiao Q, Liu Z, Wang W, Lai CH, Yang W, Yue P, Ye Q, Xiao J . TAK242 suppresses the TLR4 signaling pathway and ameliorates DCD liver IRI in rats. *Mol Med Rep*. 2019 Sep 20(3):2101-2110.
15. Bei-Yao Gao, Yi-Xing Cao, Peng-Fei Fu, Ying Xing, Dan Liang, Shan Jiang, Yu-Xiao Xie, Min Li . Optogenetics stimulates nerve reorganization in the contralesional anterolateral primary motor cortex in a mouse model of ischemic stroke *Neural Regen Res*. 2022 Jul;17(7):1535-1544.
16. Jiahong Yang, Xiaolin Sun, Ning Huang, Peng Li, Jiaqi He, Lan Jiang, Xuemei Zhang, Shu Han, Hong Xin . Entacapone alleviates acute kidney injury by inhibiting ferroptosis *FASEB J*. 2022 Jul;36(7):e22399.
17. Chi Li, Yang-Jie Ru, Quan-Yue Lin, Guang-Chun Gao, Yu-Die Yang, Xiao-Qin Zhang, Jin-Lai Gao, Shi-Hui Liu, Chu-Wei Zheng, Lin Wang, Ya-Xin Zheng, Ji-Ming Wu . Schisantherin D from *Schisandra chinensis* (Turcz.) Baill. exhibits anti-liver fibrosis capacity via modulating ETBR involved signaling, an in vitro and in vivo study *Fitoterapia*. 2022 Oct;162:105290.
18. Fengxia Liu, Mujun Li . BNIP3-mediated autophagy via the mTOR/ULK1 pathway induces primordial follicle loss after ovarian tissue transplantation *J Assist Reprod Genet*. 2023 Mar;40(3):491-508.
19. Yuanyuan Hao, Tao Song, Mingye Wang, Tongtong Li, Chi Zhao, Ting Li, Yunlong Hou, Hongjiang He . Dual targets of lethal apoptosis and protective autophagy in liver cancer with periplocymarin elicit a limited therapeutic effect *Int J Oncol*. 2023 Mar;62(3):44..
20. Mengzhi Li, Peng Wang, Jingting Li, Fei Zhou, Shixin Huang, Shaohai Qi, Bin Shu . NRP1 transduces mechanical stress inhibition via LATS1/YAP in hypertrophic scars *Cell Death Discov*. 2023 Sep 13;9(1):341..
21. Jia Li, Zheng Hu, Jiwei Zhu, Xin Lin, Xu Gao, Guixiang Lv . Antitumor Effects of Pegylated Zinc Protoporphyrin-Mediated Sonodynamic Therapy in Ovarian Cancer *Pharmaceutics*. 2023 Sep 4;15(9):2275.

Version 2025.03.12