

Protein A+G Agarose

产品编号	产品名称	包装
P2012	Protein A+G Agarose	2ml

产品简介:

- 本Protein A+G Agarose为进口分装，主要用于免疫沉淀(Immunoprecipitation, IP)或免疫共沉淀(Co-IP)，也可以用于抗体的纯化。
- Protein A+G Agarose适合于免疫沉淀所有Protein A Agarose和Protein G Agarose单独可以免疫沉淀的抗体，包括mouse IgG₁, IgG_{2a}, IgG_{2b}, IgG₃, IgA, rat IgG₁, IgG_{2a}, IgG_{2b}, IgG_{2c}, rabbit IgG, rabbit and goat polyclonal Abs, 以及human IgG₁, IgG₂, IgG₃和IgG₄。
- Protein A和Protein G都共价交联到4% agarose beads上，2ml Protein A+G Agarose中共含有约2mg重组的Protein A+G。2 ml Protein A+G Agarose共可以结合约14mg human IgG。
- Protein A+G Agarose配制在TBS溶液中，2ml中共含有0.5ml Agarose beads。
- 本Protein A+G Agarose如果用于常规的免疫沉淀，可以免疫沉淀100次。

包装清单:

产品编号	产品名称	包装
P2012	Protein A+G Agarose	1ml/管, 共2管
—	说明书	1份

保存条件:

4℃保存，一年有效。

注意事项:

- 请勿冷冻保存本产品。
- Protein A+G Agarose使用前一定要充分重悬，即充分颠倒若干次使混合均匀。
- 从蛋白样品收集开始，所有步骤中蛋白样品都必须在4℃或冰上操作。
- 为了您的安全和健康，请穿实验服并戴一次性手套操作。

使用说明:

1. 免疫沉淀(Immunoprecipitation, IP):

A. 蛋白样品的准备:

- A1. 对于10厘米细胞培养皿中的贴壁细胞，吸除细胞培养液，PBS洗涤一次，然后加入500微升至2毫升细胞裂解液裂解细胞。可以使用碧云天生产的Western及IP细胞裂解液(P0013)或各种RIPA裂解液(P0013B、P0013C、P0013D或P0013E)等进行细胞的裂解。
 - A2. 对于组织样品参考贴壁细胞使用裂解液的比例进行裂解。
 - A3. 对于悬浮细胞，离心收集细胞后，PBS洗涤一次，然后参考贴壁细胞的裂解方法进行裂解。
- 注：详细的裂解方法参考不同裂解液的详细使用方法。对于不同的培养器材，参考10厘米培养皿的裂解液的用量进行裂解。如果裂解获得的蛋白样品浓度过高，可以用裂解液或PBS适当稀释，如果蛋白样品浓度过低，在以后的裂解过程中宜适当减少裂解液的用量。

B. 去除非特异性结合(可选做):

- B1. 取200微升至1毫升蛋白样品，蛋白量约为200微克至1毫克，加入约1微克和免疫沉淀时使用的IgG种属相同的普通IgG和20微升充分重悬的Protein A+G Agarose，4℃缓慢摇动30分钟至2小时。
 - B2. 2500rpm(约1000g)离心5分钟，取上清用于后续的免疫沉淀。
- 注：所谓种属相同的IgG是指，例如后续免疫沉淀时用的是小鼠IgG，则在本步骤中可以加入normal mouse IgG，如无normal IgG可以加入其它不影响后续检测的其它mouse IgG类型的抗体。通过和normal IgG和Protein A+G Agarose的孵育，可以充分降低非特异性的结合，降低背景。

C. 免疫沉淀:

- C1. 加入0.2-2微克用于免疫沉淀的一抗，4℃缓慢摇动过夜。
- C2. 再加入20微升充分重悬的Protein A+G Agarose，4℃缓慢摇动1-3个小时。(为方便后续的洗涤操作可以把加入充分重悬的Protein A+G Agarose的量调整为40微升。)

- C3. 2500rpm(约1000g)离心5分钟, 或瞬时高速离心, 小心吸除上清, 注意宁可留下少量上清也不能吸掉Protein A+G Agarose。
- C4. 用准备蛋白样品时的裂解液或PBS洗涤沉淀5次, 裂解液或PBS的用量每次为0.5-1毫升。洗涤时离心条件和吸除上清的要求同上面的步骤C3。
- C5. 完成最后一次洗涤后, 去除上清, 加入20-40微升1XSDS-PAGE电泳上样缓冲液Vortex重悬沉淀, 瞬时高速离心把样品离心至管底。
- C6. 100℃或沸水浴处理3-5分钟, 取部分或全部样品用于SDS-PAGE电泳, 暂时不用的样品可以-20℃保存。

2. 免疫共沉淀:

参考免疫沉淀的方法进行, 但免疫共沉淀(co-IP)通常必须使用未经冻存的新鲜蛋白样品。普通的免疫沉淀虽然可以使用冻存的蛋白样品, 但也宜用新鲜的蛋白样品为佳。

使用本产品的文献:

1. Liu ZJ, Gao X, Cai Y, Yang X, Fu XL, Chen J, Zhang XB, Xin HM.
Construction of a Full-length iASPP Expression Plasmid pcDNA3.1(+)/iASPP and its biological activity.
Plasmid. 2009 Feb 2.
2. You QH, Sun GY, Wang N, Shen JL, Wang Y.
Interleukin-17F-induced pulmonary microvascular endothelial monolayer hyperpermeability via the protein kinase C pathway.
J Surg Res. 2010;162(1):110-21. Epub 2009 Feb 8.
3. Li Z, Shi K, Guan L, Cao T, Jiang Q, Yang Y, Xu C.
ROS leads to MnSOD upregulation through ERK2 translocation and p53 activation in selenite-induced apoptosis of NB4 cells.
FEBS Lett. 2010;584(11):2291-7. Epub 2010 Mar 28.
4. Guo XF, Zhu XF, Shang Y, Zhang SH, Zhen YS.
A bispecific enediyne-energized fusion protein containing ligand-based and antibody-based oligopeptides against epidermal growth factor receptor and human epidermal growth factor receptor 2 shows potent antitumor activity.
Clin Cancer Res. 2010 Apr 1;16(7):2085-94.
5. Wang D, Fang L, Li P, Sun L, Fan J, Zhang Q, Luo R, Liu X, Li K, Chen H, Chen Z, Xiao S.
The leader proteinase of foot-and-mouth disease virus negatively regulates the type I interferon pathway by acting as a viral deubiquitinase.
J Virol. 2011 Apr;85(8):3758-66.
6. Chen X, Wu X, Zhao Y, Wang G, Feng J, Li Q, Qian G.
A novel binding protein of single immunoglobulin IL-1 receptor-related molecule: Paralemmin-3.
Biochem Biophys Res Commun. 2011 Jan 28;404(4):1029-33.
7. Wang L, Lin F, Wang J, Wu J, Han R, Zhu L, Difiglia M, Qin Z.
Expression of mutant N-terminal huntingtin fragment (htt552-100Q) in astrocytes suppresses the secretion of BDNF.
Brain Res. 2012 Apr 17;1449:69-82.
8. Zhang K, Lu Y, Yang P, Li C, Sun H, Tao D, Liu Y, Zhang S, Ma Y.
HILI Inhibits TGF- β Signaling by Interacting with Hsp90 and Promoting T β R Degradation.
PLoS One. 2012;7(7):e41973. Epub 2012 Jul 27.
9. Li Y, He J, Sui S, Hu X, Zhao Y, Li N.
Clenbuterol upregulates histone demethylase JHDM2a via the β (2)-adrenoceptor/cAMP/PKA/p-CREB signaling pathway.
Cell Signal. 2012 Jul 20.
10. Ju D, He J, Zhao L, Zheng X, Yang G.
Estrogen related receptor α -induced adipogenesis is PGC-1 β -dependent.
Mol Biol Rep. 2012 Mar;39(3):3343-54.
11. Cai Y, Qiu S, Gao X, Gu SZ, Liu ZJ.
iASPP inhibits p53-independent apoptosis by inhibiting transcriptional activity of p63/p73 on promoters of proapoptotic genes.
Apoptosis. 2012 Aug;17(8):777-83.
12. Guo L, Wu WJ, Liu LD, Wang LC, Zhang Y, Wu LQ, Guan Y, Li QH.
Herpes simplex virus 1 ICP22 inhibits the transcription of viral gene promoters by binding to and blocking the recruitment of P-TEFb.
PLoS One. 2012;7(9):e45749. doi: 10.1371/journal.pone.0045749. Epub 2012 Sep 24.
13. Liu D, Zhou P, Zhang L, Gong W, Huang G, Zheng Y, He F.
HDAC1/DNMT3A-containing complex is associated with suppression of Oct4 in cervical cancer cells.
Biochemistry (Mosc). 2012 Aug;77(8):934-40. doi: 10.1134/S0006297912080159.
14. Zhao W, He H, Ren K, Li B, Zhang H, Lin Y, Shao RG.
MR-1 blocks the megakaryocytic differentiation and transition of CML from chronic phase to blast crisis through MEK dephosphorylation.
Blood Cancer J. 2013 Mar 29;3:e107. doi: 10.1038/bcj.2013.5.
15. Li H, Liu N, Wang S, Wang L, Zhao J, Su L, Zhang Y, Zhang S, Xu Z, Zhao B, Miao J.
Identification of a small molecule targeting annexin A7.
Biochim Biophys Acta. 2013 Sep;1832(9):2092-9. doi: 10.1016/j.bbamcr.2013.04.015. Epub 2013 May 4.
16. Han HY, Zhu SH, Jiang LL, Li Y, Dong H, Zhao QP, Kong CL, Huang B.
Molecular characterization and analysis of a novel calcium-dependent protein kinase from Eimeria tenella.
Parasitology. 2013 May;140(6):746-55. doi: 10.1017/S0031182012002107. Epub 2013 Feb 1.

17. Yang M, Qin Z, Zhu Y, Li Y, Qin Y, Jing Y, Liu S.
Vitamin D-binding protein in cerebrospinal fluid is associated with multiple sclerosis progression.
Mol Neurobiol. 2013 Jun;47(3):946-56. doi: 10.1007/s12035-012-8387-1. Epub 2013 Jan 22.
18. Dong QZ, Wang Y, Tang ZP, Fu L, Li QC, Wang ED, Wang EH.
Derlin-1 is overexpressed in non-small cell lung cancer and promotes cancer cell invasion via EGFR-ERK-mediated up-regulation of MMP-2 and MMP-9.
Am J Pathol. 2013 Mar;182(3):954-64. doi: 10.1016/j.ajpath.2012.11.019. Epub 2013 Jan 7.
19. Luo L, Shen L, Sun F, Ma Z.
Immunoprecipitation coupled with HPLC-MS/MS to discover the aromatase ligands from *Glycyrrhiza uralensis*.
Food Chem. 2013 May 1;138(1):315-20. doi: 10.1016/j.foodchem.2012.10.043. Epub 2012 Nov 8.
20. Liu J, Zhang X, Liu A, Liu S, Zhang L, Wu B, Hu Q.
Berberine induces apoptosis in p53-null leukemia cells by down-regulating XIAP at the post-transcriptional level.
Cell Physiol Biochem. 2013;32(5):1213-24. doi: 10.1159/000354520. Epub 2013 Nov 14.
21. Li Z, Shi K, Guan L, Jiang Q, Yang Y, Xu C.
Activation of p53 by sodium selenite switched human leukemia NB4 cells from autophagy to apoptosis.
Oncol Res. 2013;21(6):325-31. doi: 10.3727/096504014X14024160459087.
22. Li H, Huang S, Wang S, Wang L, Qi L, Zhang Y, Zhang S, Zhao B, Miao J.
Relationship between annexin A7 and integrin $\beta 4$ in autophagy.
Int J Biochem Cell Biol. 2013 Nov;45(11):2605-11. doi: 10.1016/j.biocel.2013.08.017. Epub 2013 Sep 2.
23. Xu B, Jin CH, Deng Y, Liu W, Yang TY, Feng S, Xu ZF.
Alpha-Synuclein Oligomerization in Manganese-Induced Nerve Cell Injury in Brain Slices: A Role of NO-Mediated S-Nitrosylation of Protein Disulfide Isomerase.
Mol Neurobiol. 2014 Dec;50(3):1098-110. doi: 10.1007/s12035-014-8711-z. Epub 2014 Apr 29.
24. Pan C, Fang D, Xu G, Liang J, Zhang G, Wang H, Xie L, Zhang R.
A novel acidic matrix protein, PfN44, stabilizes magnesium calcite to inhibit the crystallization of aragonite.
J Biol Chem. 2014 Jan 31;289(5):2776-87. doi: 10.1074/jbc.M113.504027. Epub 2013 Dec 3.
25. Xu YB, Yang L, Wang GF, Tong XK, Wang YJ, Yu Y, Jing JF, Feng CL, He PL, Lu W, Tang W, Zuo JP.
Benzimidazole derivative, BM601, a novel inhibitor of hepatitis B virus and HBsAg secretion.
Antiviral Res. 2014 Jul;107:6-15. doi: 10.1016/j.antiviral.2014.04.002. Epub 2014 Apr 15.
26. Zhang Z, Wu L, Wang J, Li G, Feng D, Zhang B, Li L, Yang J, Ma L, Qin H.
Opposing effects of PI3K/Akt and Smad-dependent signaling pathways in NAG-1-induced glioblastoma cell apoptosis.
PLoS One. 2014 Apr 23;9(4):e96283. doi: 10.1371/journal.pone.0096283. eCollection 2014.
27. Wang G, Chen Z, Zhang F, Jing H, Xu W, Ning S, Li Z, Liu K, Yao J, Tian X.
Blockade of PKC β protects against remote organ injury induced by intestinal ischemia and reperfusion via a p66shc-mediated mitochondrial apoptotic pathway.
Apoptosis. 2014 Sep;19(9):1342-53. doi: 10.1007/s10495-014-1008-x.
28. Zhang L, Niu Y, Zhu L, Fang J, Wang X, Wang L, Wang CC.
Different Interaction Modes for Protein-disulfide Isomerase (PDI) as an Efficient Regulator and a Specific Substrate of Endoplasmic Reticulum Oxidoreductin-1 α (Ero1 α).
J Biol Chem. 2014 Nov 7;289(45):31188-99. doi: 10.1074/jbc.M114.602961. Epub 2014 Sep 25.
29. Dong H, Wang Y, Han H, Li T, Zhao Q, Zhu S, Li L, Wu Y, Huang B.
Identification and characterization of an Eimeria-conserved protein in Eimeria tenella.
Parasitol Res. 2014 Feb;113(2):735-45. doi: 10.1007/s00436-013-3703-x. Epub 2013 Nov 24.
30. Chen Y, Hu W, Lu Y, Jiang S, Li C, Chen J, Tao D, Liu Y, Yang Y, Ma Y.
A TALEN-based specific transcript knock-down of PIWIL2 suppresses cell growth in HepG2 tumor cell.
Cell Prolif. 2014 Oct;47(5):448-56. doi: 10.1111/cpr.12120. Epub 2014 Jul 15.
31. Qiu M, Bao W, Wang J, Yang T, He X, Liao Y, Wan X.
FOXA1 promotes tumor cell proliferation through AR involving the Notch pathway in endometrial cancer.
BMC Cancer. 2014 Feb 11;14:78. doi: 10.1186/1471-2407-14-78.
32. Dong ZQ, Zhang J, Chen XM, He Q, Cao MY, Wang L, Li HQ, Xiao WF, Pan CX, Lu C, Pan MH.
Bombyx mori nucleopolyhedrovirus ORF79 is a per os infectivity factor associated with the PIF complex.
Virus Res. 2014 May 12;184:62-70. doi: 10.1016/j.virusres.2014.02.009. Epub 2014 Feb 26.
33. Li JX, Feng JM, Wang Y, Li XH, Chen XX, Su Y, Shen YY, Chen Y, Xiong B, Yang CH, Ding J, Miao ZH.
The B-Raf(V600E) inhibitor dabrafenib selectively inhibits RIP3 and alleviates acetaminophen-induced liver injury.
Cell Death Dis. 2014 Jun 5;5:e1278. doi: 10.1038/cddis.2014.241.
34. Li M, Lu Y, Hu Y, Zhai X, Xu W, Jing H, Tian X, Lin Y, Gao D, Yao J.
Salvianolic acid B protects against acute ethanol-induced liver injury through SIRT1-mediated deacetylation of p53 in rats.
Toxicol Lett. 2014 Jul 15;228(2):67-74. doi: 10.1016/j.toxlet.2014.04.011. Epub 2014 Apr 21.
35. Du Y, Bi J, Liu J, Liu X, Wu X, Jiang P, Yoo D, Zhang Y, Wu J, Wan R, Zhao X, Guo L, Sun W, Cong X, Chen L, Wang J.
3Cpro of foot-and-mouth disease virus antagonizes the interferon signaling pathway by blocking STAT1/STAT2 nuclear translocation.
J Virol. 2014 May;88(9):4908-20. doi: 10.1128/JVI.03668-13. Epub 2014 Feb 19.
36. Zhang F, Li ZL, Xu XM, Hu Y, Yao JH, Xu W, Jing HR, Wang S, Ning SL, Tian XF.
Protective effects of icariin mediated SIRT1/FOXO3 signaling pathway on intestinal ischemia/reperfusion induced acute lung injury.
Mol Med Rep. 2015 Jan;11(1):269-76. doi: 10.3892/mmr.2014.2679. Epub 2014 Oct 16.