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- Xiao-Long W, Chuan-Ping Y, Kai X, Ou-Jv Q. Selenoprotein W depletion in vitro might indicate that its main function is not as an antioxidative enzyme. *BIOCHEMISTRY-MOSCOW+* 2010;75(2):201-7.
- Wang J, Sun P, Bao Y, Liu J, An L. Cytotoxicity of single-walled carbon nanotubes on PC12 cells. *Toxicol In Vitro* 2011 Feb;25(1):242-50.
- Chen J, Xiao S, Deng Y, Du X, Yu Z. Cloning of a novel glutathione S-transferase 3 (GST3) gene and expression analysis in pearl oyster, *Pinctada martensii*. *FISH SHELLFISH IMMUN* 2011 Dec;31(6):823-30.
- Wang J, Sun P, Bao Y, Dou B, Song D, Li Y. Vitamin E renders protection to PC12 cells against oxidative damage and apoptosis induced by single-walled carbon nanotubes. *Toxicol In Vitro* 2012 Feb;26(1):32-41.
- Qian Y, Guan T, Huang M, Cao L, Li Y, Cheng H, Jin H, Yu D. Neuroprotection by the soy isoflavone, genistein, via inhibition of mitochondria-dependent apoptosis pathways and reactive oxygen induced-NF- κ B activation in a cerebral ischemia mouse model. *Neurochem Int* 2012 Jun;60(8):759-67.
- Shi H, Wang Y, Cheng Z, Ye T, Chan Z. Analysis of natural variation in bermudagrass (*Cynodon dactylon*) reveals physiological responses underlying drought tolerance. *PLoS One* 2012;7(12):e53422.
- Zhao J, Xu S, Song F, Nian L, Zhou X, Wang S. 2,3,5,4'-tetrahydroxystilbene-2-O- β -D-glucoside protects human umbilical vein endothelial cells against lysophosphatidylcholine-induced apoptosis by upregulating superoxide dismutase and glutathione peroxidase. *IUBMB Life* 2014 Oct;66(10):711-22.
- Li C, Shi L, Chen D, Ren A, Gao T, Zhao M. Functional analysis of the role of glutathione peroxidase (GPx) in the ROS signaling pathway, hyphal branching and the regulation of ganoderic acid biosynthesis in *Ganoderma lucidum*. *Fungal Genet Biol* 2015 Sep;82:168-80.
- Liu Q, Xu C, Ji G, Liu H, Mo Y, Tollerud DJ, Gu A, Zhang Q. Sublethal effects of zinc oxide nanoparticles on male reproductive cells. *Toxicol In Vitro* 2016 Sep;35:131-8.
- Sun Y, Xiu C, Liu W, Tao Y, Wang J, Qu YI. Grape seed proanthocyanidin extract protects the retina against early diabetic injury by activating the Nrf2 pathway. *Exp Ther Med* 2016 Apr;11(4):1253-1258.
- Yao C, Li G, Qian Y, Cai M, Yin H, Xiao L, Tang W, Guo F, Shi B. Protection of Pentoxifylline against Testis Injury Induced by Intermittent Hypobaric Hypoxia. *Oxid Med Cell Longev* 2016;2016:3406802.
- Zhai X, Zhang C, Zhao G, Stoll S, Ren F, Leng X. Antioxidant capacities of the selenium nanoparticles stabilized by chitosan. *J NANOBIOECHANOL* 2017 Jan 5;15(1):4.
- He JH, Yan M, Zuo H, Niu S, Yuan J, Weng SP, He J, Xu X. High reduced/oxidized glutathione ratio in infectious spleen and kidney necrosis virus-infected cells contributes to degradation of VP08R multimers. *Vet Microbiol* 2017 Aug;207:19-24.
- Zhu Z, Fan X, Lv Y, Lin Y, Wu, Zeng W. Glutamine protects rabbit spermatozoa against oxidative stress via glutathione synthesis during cryopreservation. *REPROD FERT DEVELOP* 2017 Oct;29(11):2183-2194.
- Zhu Z, Ren Z, Fan X, Pan Y, Lv S, Pan C, Lei A, Zeng W. Cysteine protects rabbit spermatozoa against reactive oxygen species-induced damages. *PLoS One* 2017 Jul 10;12(7):e0181110.
- Guo J, Xu B, Han Q, Zhou H, Xia Y, Gong C, Dai X, Li Z, Wu G. Ferroptosis: A Novel Anti-Tumor Action for Cisplatin. *Cancer Res Treat* 2017 May 10.
- Han X, Xue X, Zhao Y, Li Y, Liu W, Zhang J, Fan S. Rutin-Enriched Extract from *Coriandrum sativum* L. Ameliorates Ionizing Radiation-Induced Hematopoietic Injury. *Int J Mol Sci* 2017 Apr 29;18(5). pii: E942.
- Zhao L, Qi L, Li C, Li L, Jin L, Yuan J. SVCV impairs mitochondria complex III resulting in accumulation of hydrogen peroxide. *FISH SHELLFISH IMMUN* 2018 Apr;75:58-65.
- Lu MC, Jiao Q, Liu T, Tan SJ, Zhou HS, You QD, Jiang ZY. Discovery of a head-to-tail cyclic peptide as the Keap1-Nrf2 protein-protein interaction inhibitor with high cell potency. *Eur J Med Chem* 2018 Jan 1;143:1578-1589.
- Zhang X, Högy P, Wu X, Schmid I, Wang X, Schulze WX, Jiang D, Fangmeier A. Physiological and Proteomic Evidence for the Interactive Effects of Post-Anthesis Heat Stress and Elevated CO₂ on Wheat. *Proteomics* 2018 Dec;18(23):e1800262.
- Xu S, Zhao X, Liu F, Cao Y, Wang B, Wang X, Yin M, Wang Q, Feng X. Crucial role of oxidative stress in bactericidal effect of parthenolide against *Xanthomonas oryzae* pv. *oryzae*. *Pest Manag Sci* 2018 Dec;74(12):2716-2723.

22. Wu Z, Wang Y, Meng X, Wang X, Li Z, Qian S, Wei Y, Shu L, Ding Y, Wang P, Peng Y Total C-21 steroidal glycosides, isolated from the root tuber of *Cynanchum auriculatum* Royle ex Wight, attenuate hydrogen peroxide-induced oxidative injury and inflammation in L02 cells. *Int J Mol Med* 2018 Dec;42(6):3157-3170.
23. Fu Q, Li C, Yu L Gambogic acid inhibits spinal cord injury and inflammation through suppressing the p38 and Akt signaling pathways. *Mol Med Rep* 2018 Jan;17(1):2026-2032.
24. Dai J, Gu L, Su Y, Wang Q, Zhao Y, Chen X, Deng H, Li W, Wang G, Li K Inhibition of curcumin on influenza A virus infection and influenzal pneumonia via oxidative stress, TLR2/4, p38/JNK MAPK and NF- κ B pathways. *Int Immunopharmacol* 2018 Jan;54:177-187.
25. Chen M, Zheng J, Liu G, Xu E, Wang J, Fuqua BK, Vulpe CD, Anderson GJ, Chen H Ceruloplasmin and hephaestin jointly protect the exocrine pancreas against oxidative damage by facilitating iron efflux. *Redox Biol* 2018 Jul;17:432-439.
26. Zhang Y, Li Q, Fang M, Ma Y, Liu N, Yan X, Zhou J, Li F The Kidney Injury Induced by Short-Term PM2.5 Exposure and the Prophylactic Treatment of Essential Oils in BALB/c Mice. *Oxid Med Cell Longev* 2018 Jul 29;2018:9098627.
27. Qiao H, Ren H, Du H, Zhang M, Xiong X, Lv R Liraglutide repairs the infarcted heart: The role of the SIRT1/Parkin/mitophagy pathway. *Mol Med Rep* 2018 Mar;17(3):3722-3734.
28. Zhang QL, Wang W, Jiang Y, A-Tuya, Dongmei, Li LL, Lu ZJ, Chang H, Zhang TZ GRGM-13 comprising 13 plant and animal products, inhibited oxidative stress induced apoptosis in retinal ganglion cells by inhibiting P2RX7/p38 MAPK signaling pathway. *Biomed Pharmacother* 2018 May;101:494-500.
29. Zhang D, Li Y, Zhang T, Liu J, Jahejo AR, Yang L, Chen P, Ning G, Huo N, Ma H, Yan F, Tian W Protective effects of zinc and N-acetyl-L-cysteine supplementation against cadmium induced erythrocyte cytotoxicity in Arbor Acres broiler chickens (*Gallus gallus domesticus*). *ECOTOX ENVIRON SAFE* 2018 Nov 15;163:331-339.
30. Ding X, Sun W, Chen J IL-2 augments the sorafenib-induced apoptosis in liver cancer by promoting mitochondrial fission and activating the JNK/TAZ pathway. *Cancer Cell Int* 2018 Nov 9;18:176.
31. Deng J, Liu Q, Zhang Q, Zhang C, Liu D, Fan D, Yang H Comparative study on composition, physicochemical and antioxidant characteristics of different varieties of kiwifruit seed oil in China. *Food Chem* 2018 Oct 30;264:411-418.
32. Xu E, Chen M, Zheng J, Maimaitiming Z, Zhong T, Chen H Deletion of hephaestin and ceruloplasmin induces a serious systemic iron deficiency and disrupts iron homeostasis. *BIOCHEM BIOPH RES CO* 2018 Sep 10;503(3):1905-1910.
33. Qin QJ, Cui LQ, Li P, Wang YB, Zhang XZ, Guo ML Rhynchophylline ameliorates myocardial ischemia/reperfusion injury through the modulation of mitochondrial mechanisms to mediate myocardial apoptosis. *Mol Med Rep* 19(4):2581-2590. 2019 Apr
34. Yu HF, Duan CC, Yang ZQ, Wang YS, Yue ZP, Guo B HB-EGF Ameliorates Oxidative Stress-Mediated Uterine Decidualization Damage. *Oxid Med Cell Longev* 2019:6170936. 2019 Dec 2
35. Zhang L, Li S, Wang R, Chen C, Ma W, Cai H Cytokine augments the sorafenib-induced apoptosis in Huh7 liver cancer cell by inducing mitochondrial fragmentation and activating MAPK-JNK signalling pathway. *Biomed Pharmacother* 110:213-223. 2019 Feb
36. Sun X, Li X, Jia H, Looor JJ, Bucktrout R, Xu Q, Wang Y, Shu X, Dong J, Zuo R, Yang L, Liu G, Li X Effect of heat-shock protein B7 on oxidative stress in adipocytes from preruminant calves. *J Dairy Sci* 102(6):5673-5685. 2019 Jun
37. Alboghobeish S, Mahdavinia M, Zeidoooni L, Samimi A, Oroojan AA, Alizadeh S, Dehghani MA, Ahangarpour A, Khorsandi L Efficiency of naringin against reproductive toxicity and testicular damages induced by bisphenol A in rats. *Iran J Basic Med Sci* 22(3):315-523. 2019 Mar
38. Qiu WJ, Xu MZ, Zhu XD, Ji YH MicroRNA-27a alleviates IL-1 β -induced inflammatory response and articular cartilage degradation via TLR4/NF- κ B signaling pathway in articular chondrocytes. *Int Immunopharmacol* 76:105839. 2019 Nov
39. Lu MC, Zhao J, Liu YT, Liu T, Tao MM, You QD, Jiang ZY CPUY192018, a potent inhibitor of the Keap1-Nrf2 protein-protein interaction, alleviates renal inflammation in mice by restricting oxidative stress and NF- κ B activation. *Redox Biol* 26:101266. 2019 Sep
40. Hang Yu, Zhidian Wu, Xiaozhi Wang, Chang Gao, Run Liu, Fuxin Kang, Mingming Dai Protective effects of combined treatment with mild hypothermia and edaravone against cerebral ischemia/reperfusion injury via oxidative stress and Nrf2 pathway regulation *Int J Oncol* 2020 Aug;57(2):500-508.
41. Yangyang Chen, Jichun Tang, Yinghong Zhang, Juan Du, Yuanli Wang, Hui Yu, Yanling He Astaxanthin alleviates gestational diabetes mellitus in mice through suppression of oxidative stress *Naunyn Schmiedebergs Arch Pharmacol* 2020 Dec;393(12):2517-2527.
42. Qiushi Xu, Yunhui Fan, Juan J Looor, Yusheng Liang, Xudong Sun, Hongdou Jia, Chenxu Zhao, Chuang Xu Adenosine 5'-monophosphate-activated protein kinase ameliorates bovine adipocyte oxidative stress by inducing antioxidant responses and autophagy *J Dairy Sci* 2021 Apr;104(4):4516-4528.
43. Li Ma, Chengxu Li, Shuai Lian, Bin Xu, Hongming Lv, Yanzhi Liu, Jingjing Lu, Hong Ji, Shize Li, Jingru Guo, Huanmin Yang Procyanidin B2 alleviates liver injury caused by cold stimulation through Sonic hedgehog signalling and autophagy *J Cell Mol Med* 2021 Aug;25(16):8015-8027.
44. Yan He, Yuanzhu Sun, Jun Peng *Circ* 0114428 Regulates Sepsis-Induced Kidney Injury by Targeting the miR-495-3p/CRBN Axis *Inflammation* 2021 Aug;44(4):1464-1477.
45. Kun Ma, Hao-Yu Wu, Sheng-Yuan Wang, Bai-Xiang Li The Keap1/Nrf2-ARE signaling pathway is involved in atrazine induced dopaminergic neurons degeneration via microglia activation *Ecotoxicol Environ Saf* 2021 Dec 15;226:112862.
46. Songge Deng, Yizhao Wu, Hanqi Duan, Jo-Anne E Cavanagh, Xiuhong Wang, Jiangping Qiu, Yinsheng Li Toxicity assessment of earthworm exposed to arsenate using oxidative stress and burrowing behavior responses and an integrated biomarker index *Sci Total Environ* 2021 Dec 15;800:149479.
47. Juanping Yin, Xinke Chen Edaravone prevents high glucose-induced injury in retinal Müller cells through thioredoxin1 and the PGC-1 α /NRF1/TFAM pathway *Pharm Biol* 2021 Dec;59(1):1233-1244.
48. Yike Yang, Ping Xu, Fangyu Zhu, Jiujiang Liao, Yue Wu, Mingyu Hu, Huijia Fu, Juan Qiao, Li Lin, Biao Huang, Huili Jin, Xiyao Liu, Yangxi Zheng, Li Wen, Richard Saffery, Mark D Kilby, Jianying Yan, Louise C Kenny, Hongbo Qi, Chao Tong, Philip N Baker The Potent Antioxidant MitoQ Protects Against Preeclampsia During Late Gestation but Increases the Risk of Preeclampsia When Administered in Early Pregnancy *Antioxid Redox Signal* 2021 Jan 10;34(2):118-136.
49. Xiuzhen Hou, Junfeng Zhang, Hehong Ma, Ming Li, Pei Wang Hypoxia-reoxygenation treatment attenuates gestational diabetes mellitus *Endocr Connect* 2021 Jan;10(1):84-91.
50. Runpu Li, Yingxin Qu, Xiaoqi Li, Ye Tao, Qinghua Yang, Junyi Wang, Yumei Diao, Qian Li, Yifan Fang, Yifei Huang, Liqiang Wang

- Molecular Hydrogen Attenuated N-methyl-N-Nitrosourea Induced Corneal Endothelial Injury by Upregulating Anti-Apoptotic Pathway Invest Ophthalmol Vis Sci 2021 Jul 1;62(9):2.
51. Yongshi Ma, Min Xuan, Yunqing Dong, Hong Wang, Jianfang Huang, Biao Cheng, Xiao Cui Platelet-rich plasma protects human melanocytes from oxidative stress and ameliorates melanogenesis induced by UVB irradiation Biosci Biotechnol Biochem 2021 Jun 24;85(7):1686-1696.
 52. Xiao-Huan Zhang, Chen-Cong Zhou, Chong-Yong Li, Yao Hua, Kang Li, Ping Wei, Ming-Fang He Isoliquiritin exert protective effect on telencephalon infarction injury by regulating multi-pathways in zebrafish model of ischemic stroke Phytomedicine 2021 Mar;83:153469.
 53. Xiao Cui, Yongshi Ma, Hong Wang, Jianfang Huang, Linlin Li, Jianbing Tang, Biao Cheng The Anti-photoaging Effects of Pre- and Post-treatment of Platelet-rich Plasma on UVB-damaged HaCaT Keratinocytes Photochem Photobiol 2021 May;97(3):589-599.
 54. QingJian Liang, WenNa Dong, FeiFei Wang, WeiNa Wang, Jian Zhang, Xing Liu Ficus hirta Vahl. promotes antioxidant enzyme activity under ammonia stress by inhibiting miR-2765 expression in Penaeus vannamei Ecotoxicol Environ Saf 2021 Nov 15;228:112989.
 55. Jingtao Jiang, Chengji Dong, Liang Zhai, Junsheng Lou, Jie Jin, Sheng Cheng, Zhuliu Chen, Xiaoshan Guo, Damu Lin, Jian Ding, Weiyang Gao Paeoniflorin Suppresses TBHP-Induced Oxidative Stress and Apoptosis in Human Umbilical Vein Endothelial Cells via the Nrf2/HO-1 Signaling Pathway and Improves Skin Flap Survival Front Pharmacol 2021 Nov 4;12:735530.
 56. Hui Zhu, Chunyi Lu, Fei Gao, Zhixiang Qian, Yu Yin, Shidong Kan, Daijie Chen Selenium-enriched Bifidobacterium longum DD98 attenuates irinotecan-induced intestinal and hepatic toxicity in vitro and in vivo Biomed Pharmacother 2021 Nov;143:112192.
 57. Qiancheng Mu, Yuxuan Zhang, Long Gu, Stefan T Gerner, Xiancheng Qiu, Qianke Tao, Jinwei Pang, Ghosh Dipritu, Lifang Zhang, Shigang Yin, Yong Jiang, Jianhua Peng Transcriptomic Profiling Reveals the Antiapoptosis and Antioxidant Stress Effects of Fos in Ischemic Stroke Front Neurol 2021 Oct 21;12:728984.
 58. Yuxia Feng, Yuxin Hu, Pengpeng Fang, Xiangjun Zuo, Jinxiong Wang, Jiana Li, Wei Qian, Jiaqin Mei Silicon Alleviates the Disease Severity of Sclerotinia Stem Rot in Rapeseed Front Plant Sci 2021 Sep 13;12:721436.
 59. Yaping Zhan, Minxia Zhu, Shang Liu, Jiayue Lu, Zhaohui Ni, Hong Cai, Weiming Zhang MicroRNA-93 inhibits the apoptosis and inflammatory response of tubular epithelial cells via the PTEN/AKT/mTOR pathway in acute kidney injury Mol Med Rep 2021 Sep;24(3):666.
 60. Yue Wu, Cunlin Gu, L U Huang, Yuanqing Zhao, Yanjun Tang, Hongqian Zhao, Qiong Wu Hypoxia preconditioning improves structure and function of astrocytes mitochondria via PGC-1 α /HIF signal J Biosci 2021;46:7.
 61. Guanning Huang, Jiankun Zang, Lizhen He, Huili Zhu, Jiarun Huang, Zhongwen Yuan, Tianfeng Chen, Anding Xu Bioactive Nanoenzyme Reverses Oxidative Damage and Endoplasmic Reticulum Stress in Neurons under Ischemic Stroke ACS Nano 2022 Jan 25;16(1):431-452.
 62. Bingxin Zheng, Yixin Zheng, Ningning Zhang, Yi Zhang, Baodong Zheng Rhoifolin from Plumula Nelumbinis exhibits anti-cancer effects in pancreatic cancer via AKT/JNK signaling pathways Sci Rep 2022 Apr 5;12(1):5654.
 63. Lili Song, Congjun Li, Fei Wu, Shicui Zhang Dietary intake of diosgenin delays aging of male fish Nothobranchius guentheri through modulation of multiple pathways that play prominent roles in ROS production Biogerontology 2022 Apr;23(2):201-213.
 64. Jie Song, Jinxiang Wang SIRT3 regulates bronchial epithelium apoptosis and aggravates airway inflammation in asthma Mol Med Rep 2022 Apr;25(4):144.
 65. Shaofang Zhang, Yonghui Li, Si Sun, Ling Liu, Xiaoyu Mu, Shuhu Liu, Menglu Jiao, Xinzhu Chen, Ke Chen, Huizhen Ma, Tuo Li, Xiaoyu Liu, Hao Wang, Jianning Zhang, Jiang Yang, Xiao-Dong Zhang Single-atom nanozymes catalytically surpassing naturally occurring enzymes as sustained stitching for brain trauma Nat Commun 2022 Aug 12;13(1):4744.
 66. Yinhong Niu, Jiale Liao, Haitao Zhou, Chih-Chen Wang, Lei Wang, Yanli Fan Flavonoids from Lycium barbarum Leaves Exhibit Anti-Aging Effects through the Redox-Modulation Molecules 2022 Aug 3;27(15):4952.
 67. Bingru Nan, Xia Sun, Jie Zhang, Qingyu Huang, Xi Zhang, Yanbo Li, Junchao Duan, Rui Chen, Zhiwei Sun, Heqing Shen Accumulated oxidative stress risk in HUVECs by chronic exposure to non-observable acute effect levels of PM2.5 Toxicol In Vitro 2022 Aug;82:105376.
 68. Lili Song, Kailiang Leng, Kun Xiao, Shicui Zhang Administration of krill oil extends lifespan of fish Nothobranchius guentheri via enhancement of antioxidant system and suppression of NF- κ B pathway Fish Physiol Biochem 2022 Aug;48(4):1057-1073.
 69. Dongyu Lei, Biao Li, Zuribia Isa, Xiaoqiang Ma, Baojian Zhang Hypoxia-elicited cardiac microvascular endothelial cell-derived exosomal miR-210-3p alleviate hypoxia/reoxygenation-induced myocardial cell injury through inhibiting transferrin receptor 1-mediated fe Tissue Cell 2022 Dec;79:101956.
 70. Yuqiong Li, Huiyu Wen, Yuwei Yang, Zhengwei Zhao, Haihui Gao, Hongbing Li, Meizhou Huang Potential prognostic markers of retained placenta in dairy cows identified by plasma metabolomics coupled with clinical laboratory indicators Vet Q 2022 Dec;42(1):199-212.
 71. Guang-Hui Chen, Chang-Chun Song, Tao Zhao, Christer Hogstrand, Xiao-Lei Wei, Wu-Hong Lv, Yu-Feng Song, Zhi Luo Mitochondria-Dependent Oxidative Stress Mediates ZnO Nanoparticle (ZnO NP)-Induced Mitophagy and Lipotoxicity in Freshwater Teleost Fish Environ Sci Technol 2022 Feb 15;56(4):2407-2420.
 72. Guang-Hui Chen, Chang-Chun Song, Kostas Pantopoulos, Xiao-Lei Wei, Hua Zheng, Zhi Luo Mitochondrial oxidative stress mediated Fe-induced ferroptosis via the NRF2-ARE pathway Free Radic Biol Med 2022 Feb 20:180:95-107.
 73. Binbing Liu, Yang Tian, Yuchen Li, Pei Wu, Yongzhi Zhang, Jiaolin Zheng, Huaizhang Shi ACEA Attenuates Oxidative Stress by Promoting Mitophagy via CB1R/Nrf1/PINK1 Pathway after Subarachnoid Hemorrhage in Rats Oxid Med Cell Longev 2022 Feb 24:2022:1024279.
 74. Lei Zhang, Cundong Fan, Hua-Chen Jiao, Qian Zhang, Yue-Hua Jiang, Jie Cui, Yang Liu, Yong-Hao Jiang, Juan Zhang, Meng-Qi Yang, Yan Li, Yi-Tao Xue Calycosin Alleviates Doxorubicin-Induced Cardiotoxicity and Pyroptosis by Inhibiting NLRP3 Inflammasome Activation Oxid Med Cell Longev 2022 Jan 5:2022:1733834.
 75. Zi-Run Jin, Ya-Lei Cao, Zhi-Chao Luo, Qian-Cheng Zhao, Yu Xi, Jia-Ming Weng, Zhe Zhang, Hui Jiang Therapeutic Effects of Xianlu Oral Solution on Rats with Oligoasthenozoospermia through Alleviating Apoptosis and Oxidative Stress Evid Based Complement Alternat Med 2022 Jun 18:2022:1269530.
 76. Cuicui Zhang, Hai Niu, Chengyu Wan, Xiuxian Yu, Guang Xin, Yuda

- Zhu, Zeliang Wei, Fan Li, Yilan Wang, Kun Zhang, Shiyi Li, Yuman Dong, Yangying Li, Wen Huang Drug D, a Diosgenin Derive, Inhibits L-Arginine-Induced Acute Pancreatitis through Meditating GSDMD in the Endoplasmic Reticulum via the TXNIP/HIF-1 α Pathway *Nutrients* 2022 Jun 22;14(13):2591.
77. Disi Deng, Jin Yan, Wanjing Li, Yeke Wu, Keming Wu Protective Effect of XinJiaCongRongTuSiZiWan on the Reproductive Toxicity of Female Rats Induced by Triptolide *Evid Based Complement Alternat Med* 2022 Jun 6:2022:3642349.
78. Xiliang Du, Chiara Di Malta, Zhiyuan Fang, Taiyu Shen, Xiaodi Niu, Meng Chen, Bo Jin, Hao Yu, Lin Lei, Wenwen Gao, Yuxiang Song, Zhe Wang, Chuang Xu, Zhijun Cao, Guowen Liu, Xinwei Li Nuciferine protects against high-fat diet-induced hepatic steatosis and insulin resistance via activating TFEB-mediated autophagy-lysosomal pathway *Acta Pharm Sin B* 2022 Jun;12(6):2869-2886.
79. Lili Song, Fei Wu, Congjun Li, Shicui Zhang Dietary intake of GDF11 delays the onset of several biomarkers of aging in male mice through anti-oxidant system via Smad2/3 pathway *Biogerontology* 2022 Jun;23(3):341-362.
80. Rui Wang, Jinjin Pan, Jinzhen Han, Miaomiao Gong, Liang Liu, Yunlong Zhang, Ying Liu, Dingyou Wang, Qing Tang, Na Wu, Lin Wang, Jinsong Yan, Hua Li, Yuhui Yuan Melatonin Attenuates Dasatinib-Aggravated Hypoxic Pulmonary Hypertension via Inhibiting Pulmonary Vascular Remodeling *Front Cardiovasc Med* 2022 Mar 24:9:790921.
81. Xinghui Wang, Mengyao Zhu, Juan J Looor, Qianming Jiang, Yiwei Zhu, Wei Li, Xiliang Du, Yuxiang Song, Wenwen Gao, Lin Lei, Jianguo Wang, Guowen Liu, Xinwei Li Propionate alleviates fatty acid-induced mitochondrial dysfunction, oxidative stress, and apoptosis by upregulating PPAR γ coactivator 1 α in hepatocytes *J Dairy Sci* 2022 May;105(5):4581-4592.
82. Shuqin Du, Hanqiang Shi, Lie Xiong, Ping Wang, Yanbo Shi Canagliflozin mitigates ferroptosis and improves myocardial oxidative stress in mice with diabetic cardiomyopathy *Front Endocrinol (Lausanne)* 2022 Oct 14:13:1011669.
83. Mengrong Miao, Yangyang Wang, Shuang Zeng, Yaqian Han, Ruilou Zhu, Pengfei Yu, Yitian Yang, Ningning Fu, Ningning Li, Mingyang Sun, Jiaqiang Zhang Identification and Validation of Ferroptosis-Related Genes in Sevoflurane-Induced Hippocampal Neurotoxicity *Oxid Med Cell Longev* 2022 Oct 4:2022:4435161.
84. Tianyi Wang, Tingting Zhou, Mingming Xu, Shuo Wang, Anqi Wu, Mingyang Zhang, You Lang Zhou, Jiahai Shi Platelet membrane-camouflaged nanoparticles carry microRNA inhibitor against myocardial ischaemia-reperfusion injury *J Nanobiotechnology* 2022 Oct 4;20(1):434.
85. Guang-Hui Liu, Dian-Guang Zhang, Xi-Jun Lei, Xiao-Ying Tan, Chang-Chun Song, Hua Zheng, Zhi Luo Effects of Dietary Selenium and Oxidized Fish Oils on Intestinal Lipid Metabolism and Antioxidant Responses of Yellow Catfish *Pelteobagrus fulvidraco* Antioxidants (Basel) 2022 Sep 26;11(10):1904.
86. Yangying Li, Yuda Zhu, Shiyi Li, Yuman Dong, Chengyu Wan, Xiuxian Yu, Guang Xin, Zeliang Wei, Fan Li, Yilan Wang, Kun Zhang, Qingqiu Chen, Cuicui Zhang, E Wen, Hai Niu, Wen Huang Deoxyarbutin attenuates severe acute pancreatitis via the HtrA2/PGC-1 α pathway *Free Radic Res* 2022 Sep-Oct;56(9-10):651-665.

Version 2024.03.12